



# KAISER PERMANENTE®

<b>DOCUMENT NUMBER:</b>
<b>DOCUMENT TITLE:</b>
<b>DOCUMENT NOTES:</b>

<b>LOCATION:</b>	<b>VERSION:</b>
<b>DOC TYPE:</b>	<b>STATUS:</b>

<b>EFFECTIVE DATE:</b>	<b>NEXT REVIEW DATE:</b>
<b>RELEASE DATE:</b>	<b>EXPIRATION DATE:</b>


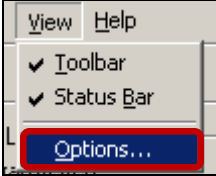
<b>AUTHOR:</b>	<b>PREVIOUS NUMBER:</b>
<b>OWNER:</b>	<b>CHANGE NUMBER:</b>

## Transferring / Tracking Specimens

**Purpose** This procedure provides instructions for transferring and tracking specimens from their current location to the testing laboratory location.

**Scope** This procedure is intended for all users.

**Adding Temperature to the Transfer List Fields** Follow the steps below to add **temperature** to the Transfer List fields.

Step	Action	Icon
1	Click on the <b>Transfer Specimen</b> icon from the Applications bar (Appbar) to launch the application.	
2	Click on <b>View</b> on the menu bar, and then select <b>Options</b> .  	
3	Check the box for <b>Filter by transfer temperature</b> .	
4	Select a default <b>Printer</b> .	
5	Click on <b>Save</b> .	

*Continued on next page*


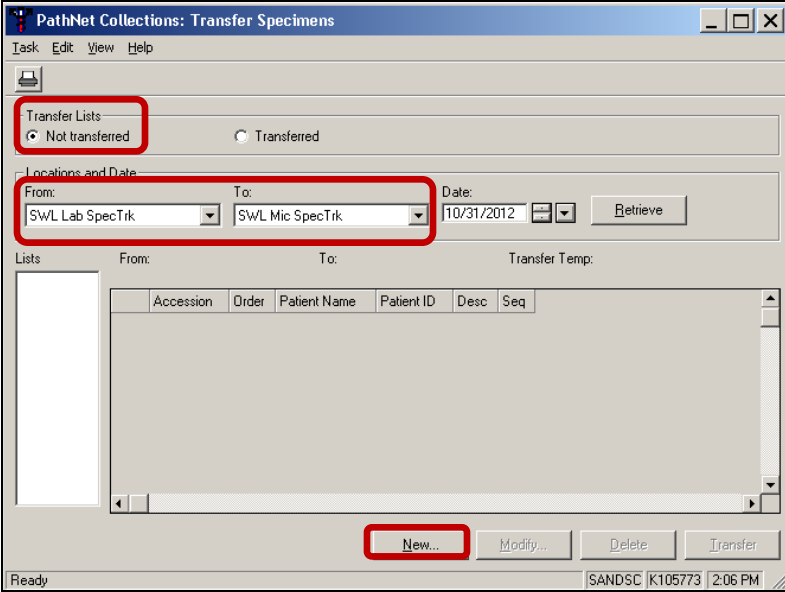
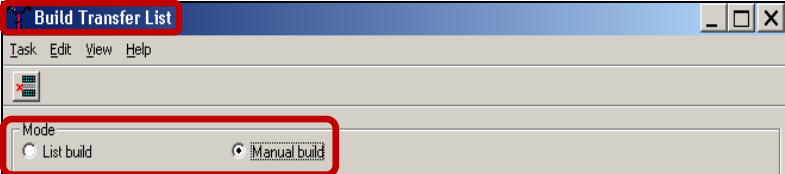
## Transferring / Tracking Specimens, Continued

### Creating a Transfer List (Manual Build mode)

Follow the steps below to create a transfer list (**Manual Build mode**).

**Note:**

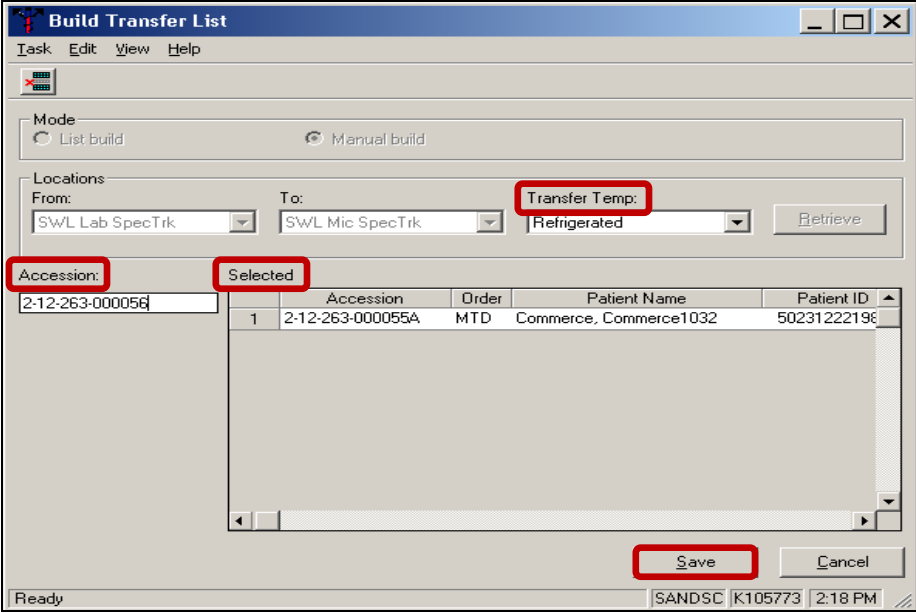
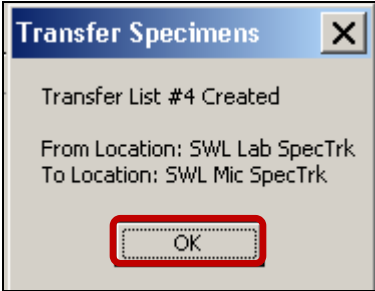
Specimens must be logged-in before they can be transferred.

Step	Action	
1	Click on the <b>Transfer Specimen icon</b> from the Appbar to launch the application.	
2	<ul style="list-style-type: none"> <li>• The <b>Transfer Specimens</b> window opens.</li> <li>• Verify that <b>Not transferred</b> is selected.</li> <li>• Using the drop-down window, select the <b>From</b> Location.</li> <li>• Using the drop-down window, select the <b>To</b> Location.                             <ul style="list-style-type: none"> <li>• The <b>To</b> Location is the <b>performing</b> location.</li> </ul> </li> <li>• Click on the <b>New</b> button at the bottom right of the screen.</li> </ul>	
3	<ul style="list-style-type: none"> <li>• The <b>Build Transfer List</b> dialog window opens.</li> <li>• Under <b>Mode</b>, select <b>Manual Build</b>.</li> </ul>	

*Continued on next page*

## Transferring / Tracking Specimens, Continued

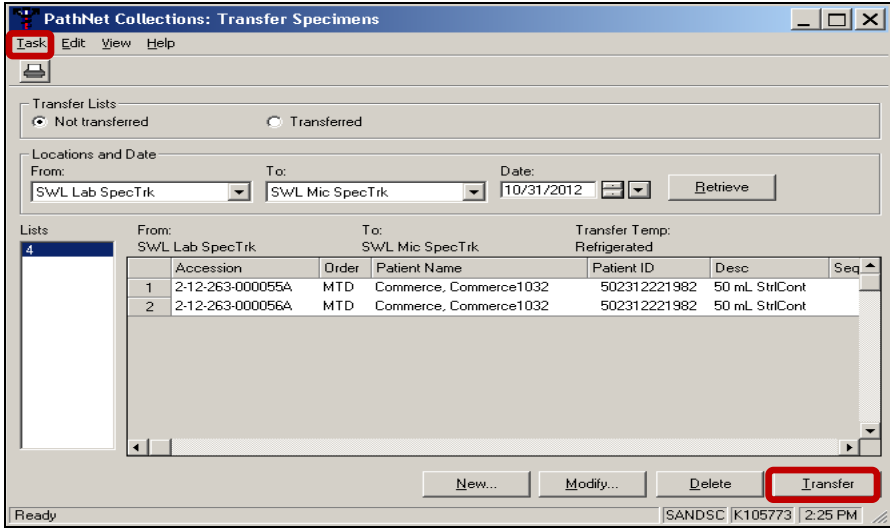
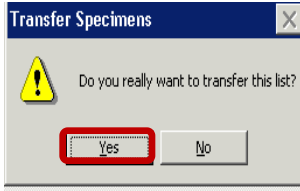
### Creating a Transfer List (Manual Build mode), continued

Step	Action
4	<ul style="list-style-type: none"> <li>• Select the <b>Transfer Temperature</b>.</li> <li>• Click in the <b>Accession</b> field.</li> <li>• Scan or type the accession number of the specimen to be added to the transfer list. Press <b>Enter</b> to move the accession number from the <b>Accession</b> window to the <b>Selected</b> window. Repeat the step for each specimen to be added to the Transfer List.</li> <li>• Click on <b>Save</b>.</li> </ul> 
5	<ul style="list-style-type: none"> <li>• The <b>Transfer Specimens</b> dialog window is displayed.</li> <li>• Click on the <b>OK</b> button to close the window.</li> </ul> 

*Continued on next page*

## Transferring / Tracking Specimens, Continued

### Creating a Transfer List (Manual Build mode), continued


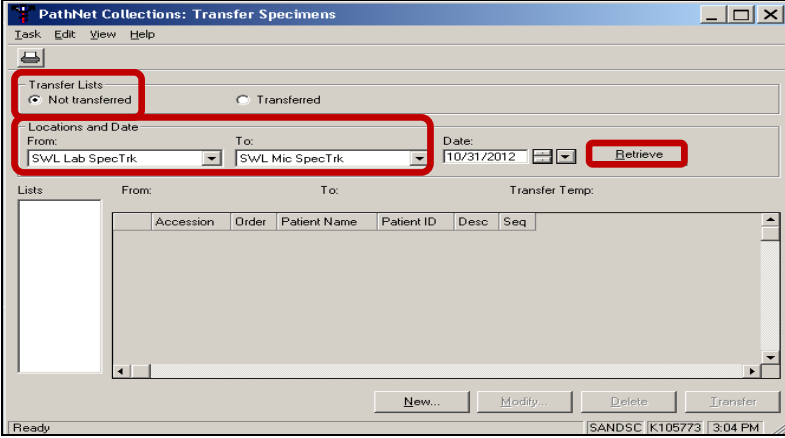
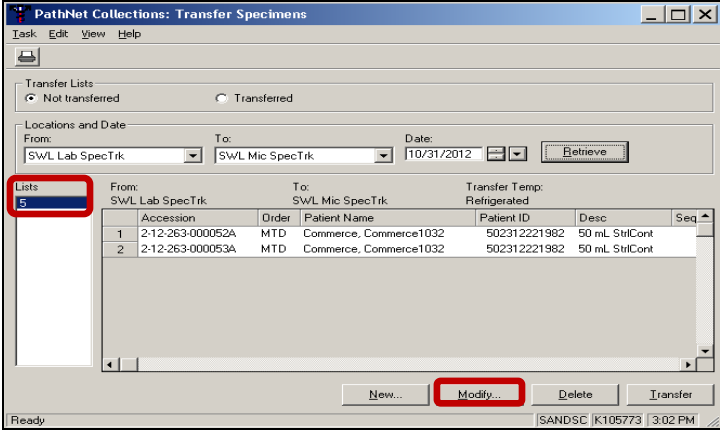
Step	Action
6	<ul style="list-style-type: none"> <li>If additional specimens will be added to the list at a later time, click on <b>T</b>ask, then on <b>E</b>xit on the Menu Bar to exit the <b>Specimen Transfer</b> window. Proceed to <b>Modifying a Transfer List</b>.</li> </ul> <p><b>OR</b></p> <ul style="list-style-type: none"> <li>To close the list so no additional specimens can be added, click on <b>T</b>ransfer.*</li> </ul> <p><i>*Note:                      Any open Transfer Lists are to be closed a few minutes prior to the established Courier pick-up time, or at the end of the business day if you are working at an MOB.</i></p> 
7	<ul style="list-style-type: none"> <li>Click on the <b>Y</b>es button to transfer the list.</li> <li>Select a printer when prompted.</li> </ul> 

Continued on next page

## Transferring / Tracking Specimens, Continued

### Modifying a Transfer List

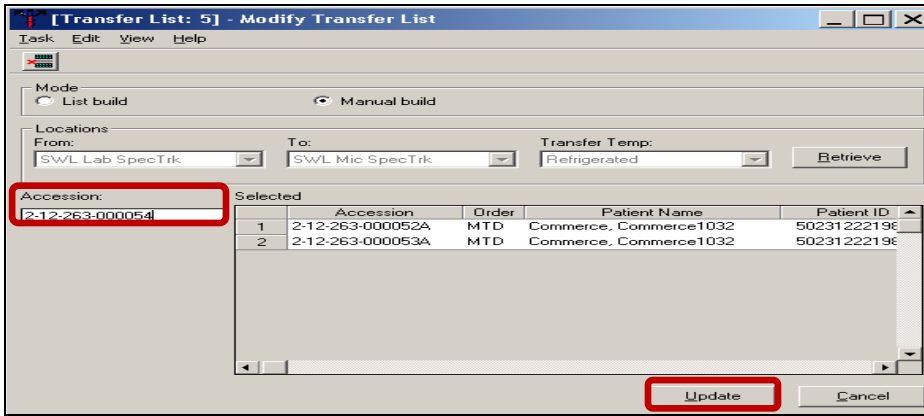
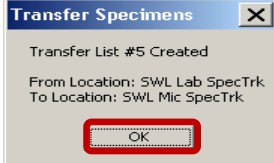
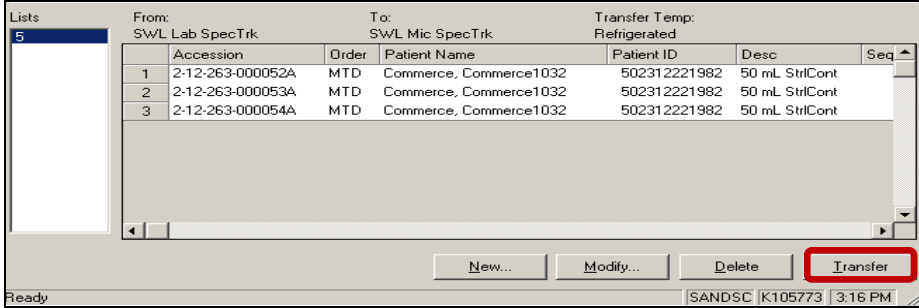
Follow the steps below to modify a transfer list.

Step	Action	Icon
1	<ul style="list-style-type: none"> <li>Click on the <b>Transfer Specimen</b> icon from the Appbar to launch the application.</li> </ul>	
2	<ul style="list-style-type: none"> <li>The <b>Transfer Specimens</b> window opens.</li> <li>Verify that <b>Not transferred</b> is selected.</li> <li>Using the drop-down window, select the <b>From</b> Location.</li> <li>Using the drop-down window, select the <b>To</b> Location.</li> <li>Verify the <b>Date</b>.</li> <li>Click on <b>Retrieve</b>.</li> </ul> 	
3	<ul style="list-style-type: none"> <li>Select the <b>Transfer List Number</b> to which specimens will be added.</li> <li>Click on the <b>Modify</b> button.</li> </ul> 	

*Continued on next page*

## Transferring / Tracking Specimens, Continued

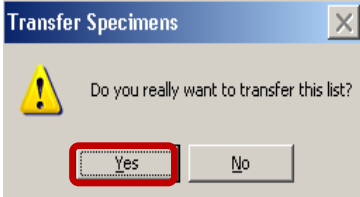
### Modifying a Transfer List, continued

Step	Action
4	<ul style="list-style-type: none"> <li>• In the Accession field, scan the <b>label barcode</b> for the specimen to be added to the selected transfer list. Press <b>Enter</b>.</li> <li>• Repeat the step for all specimens to be added to the list at this time.</li> <li>• Click on <b>Update</b> to save the additions.</li> </ul> 
5	<p>The <b>Transfer Specimen</b> window appears showing the transfer list number.                      Click on <b>OK</b>.</p> 
6	<p>Click on <b>Transfer</b> to close the list. No additional specimens can be added.</p> 

Continued on next page



## Transferring / Tracking Specimens, Continued

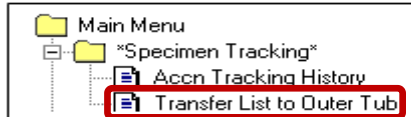
### Modifying a Transfer List, continued

Step	Action
7	<ul style="list-style-type: none"> <li>Click on the <b>Yes</b> button to transfer the list.</li> <li>Select a printer when prompted.</li> </ul> 

### Associating the Transfer List with Outer Tub

Follow the steps below to associate the Transfer List with Outer Tub Tracking Number.

Step	Action	Icon
1	Access the <b>Explorer Menu</b> icon on the Cerner AppBar.	
2	<ul style="list-style-type: none"> <li>Double-click on <b>Main Menu</b>.</li> <li>Click on the + sign to the left of the <b>Specimen Tracking</b> folder.</li> </ul> 	
3	Click on <b>Transfer List to Outer Tub</b> .	

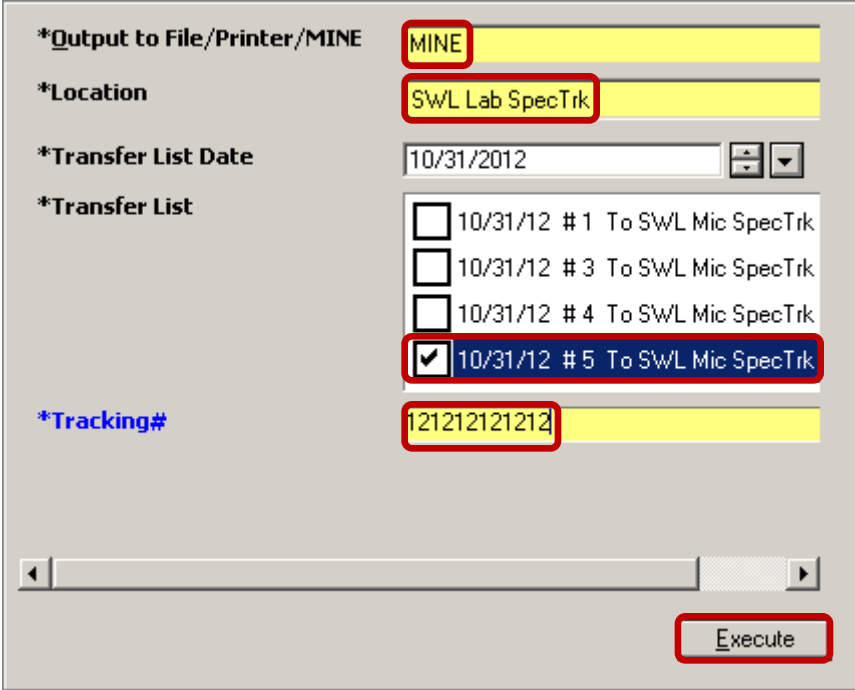


Continued on next page



## Transferring / Tracking Specimens, Continued

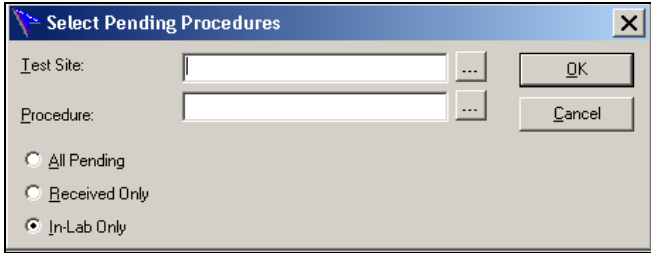

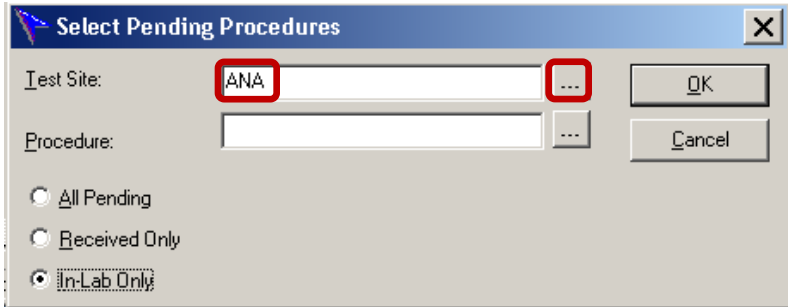
### Associating the Transfer List with Outer Tub, continued

Step	Action
4	<ul style="list-style-type: none"> <li>• Enter the printer name in the <b>*Output to file/Printer/MINE</b> field.</li> <li>• Select the <b>*Location</b> from the drop-down menu.</li> <li>• Select the <b>*Transfer List(s)</b> to be associated with the outer tub.</li> <li>• Click in the <b>*Tracking#</b> field. Scan the <b>barcode</b> for the outer tub.</li> <li>• Click on <b>Execute</b>. The <b>Transfer List to Outer Tub</b> report prints.</li> </ul> <p><b>Note:</b>                      If MINE is selected in the <b>Output...</b> field, the report will not automatically print when Execute is clicked.</p> 
5	Secure the Outer Tub in preparation for Courier pickup.

*Continued on next page*

## Transferring / Tracking Specimens, Continued

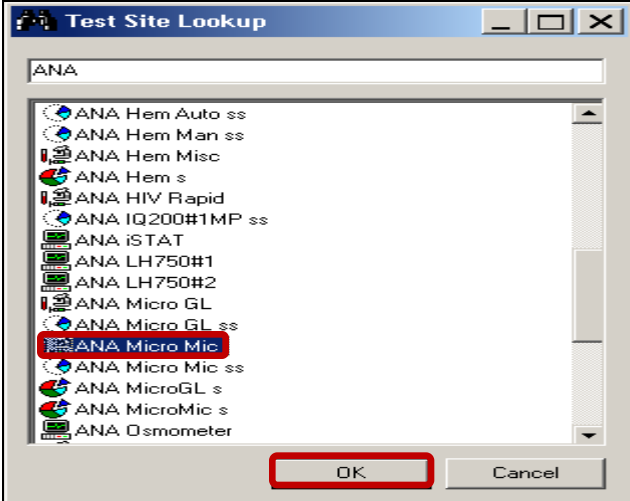
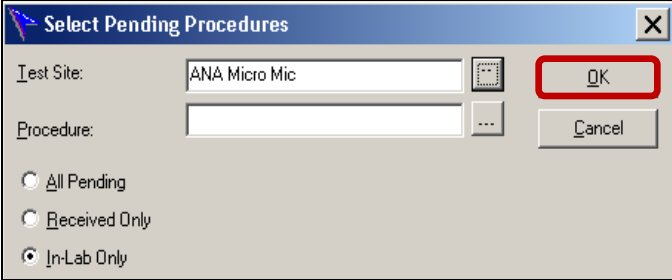
**Changing the Designated Test Site** Follow the steps below to change the designated test site for at test or group of tests.

Step	Action	Icon
1	<ul style="list-style-type: none"> <li>Click on the <b>Pending Inquiry</b> from the Appbar to launch the application.</li> <li>The <b>Select Pending Procedures</b> window opens.</li> </ul> 	
2	<ul style="list-style-type: none"> <li>Enter the <b>Test Site</b> (site abbreviation) and click on the <b>ellipses</b>.</li> <li>Enter the <b>Procedure</b> (if desired).</li> <li>Select the <b>Order Status</b> for the order(s) to be transferred.                             <ul style="list-style-type: none"> <li><b>Test Site</b> is the site abbreviation.</li> <li><b>Procedure</b> is the test name.</li> <li><b>All Pending</b> displays all pending orders for the test site and/or procedure.</li> <li>Both <b>Received Only</b> and <b>In-Lab Only</b> display all pending orders with an In-Lab status.</li> </ul> </li> </ul> 	

*Continued on next page*

## Transferring / Tracking Specimens, Continued

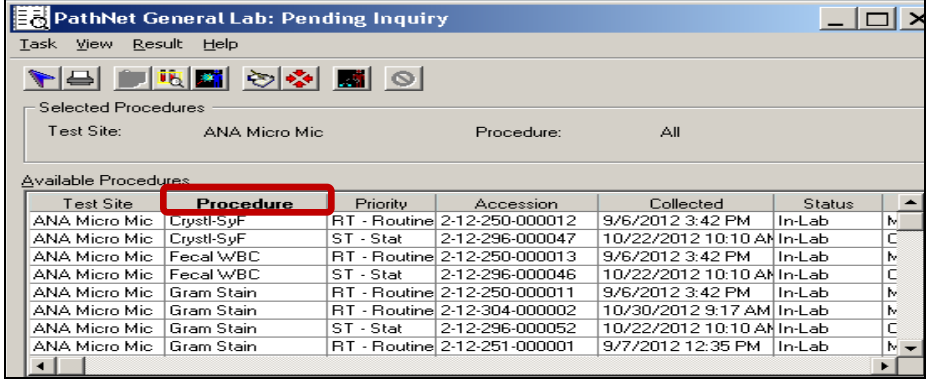
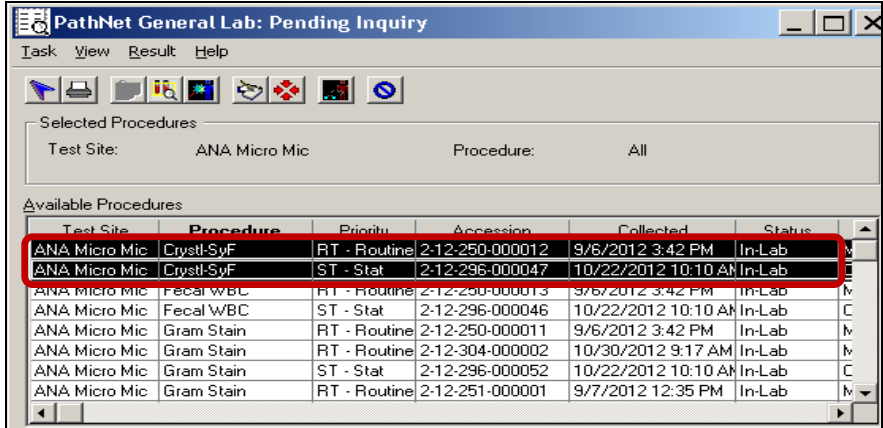
### Changing the Designated Test Site, continued

Step	Action
3	<p>The <b>Test Site Lookup</b> window opens.</p> <ul style="list-style-type: none"><li>• Select the specific site from the drop-down list.</li><li>• Click on <b>OK</b>.</li></ul>  <p>The <b>Test Site</b> box is populated with the selected specific site.</p> <ul style="list-style-type: none"><li>• Click on <b>OK</b>.</li></ul> 

*Continued on next page*

## Transferring / Tracking Specimens, Continued

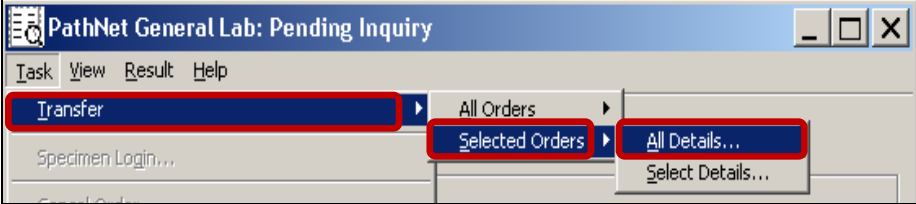

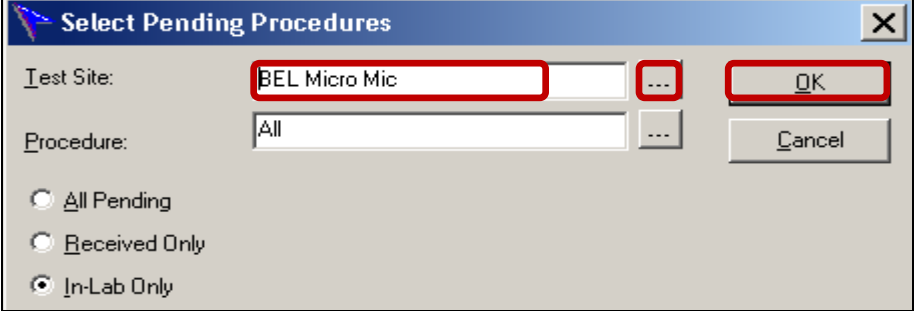
### Changing the Designated Test Site, continued

Step	Action						
4	<p>The <b>Pending Inquiry</b> window opens.</p> <ul style="list-style-type: none"> <li>Clicking on the column header will sort data in that column.</li> </ul> 						
5	<p>Select the order or orders to be transferred.</p> <ul style="list-style-type: none"> <li>For a single accession, click on the <b>row</b> containing that accession.</li> <li>For multiple orders listed in sequence:                     <table border="1" data-bbox="532 1087 1414 1398"> <thead> <tr> <th data-bbox="540 1094 760 1125">If...</th> <th data-bbox="764 1094 1406 1125">Then...</th> </tr> </thead> <tbody> <tr> <td data-bbox="540 1131 760 1262">the orders are in sequence</td> <td data-bbox="764 1131 1406 1262"> <ul style="list-style-type: none"> <li>Click on the <b>first order</b> in the sequence.</li> <li>Press and hold the <b>Shift</b> key.</li> <li>Click on the <b>last order</b> in the sequence.</li> <li>Release the <b>Shift</b> key.</li> </ul> </td> </tr> <tr> <td data-bbox="540 1268 760 1398">the orders are not in sequence</td> <td data-bbox="764 1268 1406 1398"> <ul style="list-style-type: none"> <li>Click on the <b>first order</b> to be transferred.</li> <li>Press and hold the <b>Ctrl</b> key.</li> <li>Click on any <b>additional orders</b> to be transferred.</li> <li>Release the <b>Ctrl</b> key.</li> </ul> </td> </tr> </tbody> </table> </li> </ul> 	If...	Then...	the orders are in sequence	<ul style="list-style-type: none"> <li>Click on the <b>first order</b> in the sequence.</li> <li>Press and hold the <b>Shift</b> key.</li> <li>Click on the <b>last order</b> in the sequence.</li> <li>Release the <b>Shift</b> key.</li> </ul>	the orders are not in sequence	<ul style="list-style-type: none"> <li>Click on the <b>first order</b> to be transferred.</li> <li>Press and hold the <b>Ctrl</b> key.</li> <li>Click on any <b>additional orders</b> to be transferred.</li> <li>Release the <b>Ctrl</b> key.</li> </ul>
If...	Then...						
the orders are in sequence	<ul style="list-style-type: none"> <li>Click on the <b>first order</b> in the sequence.</li> <li>Press and hold the <b>Shift</b> key.</li> <li>Click on the <b>last order</b> in the sequence.</li> <li>Release the <b>Shift</b> key.</li> </ul>						
the orders are not in sequence	<ul style="list-style-type: none"> <li>Click on the <b>first order</b> to be transferred.</li> <li>Press and hold the <b>Ctrl</b> key.</li> <li>Click on any <b>additional orders</b> to be transferred.</li> <li>Release the <b>Ctrl</b> key.</li> </ul>						

Continued on next page

## Transferring / Tracking Specimens, Continued

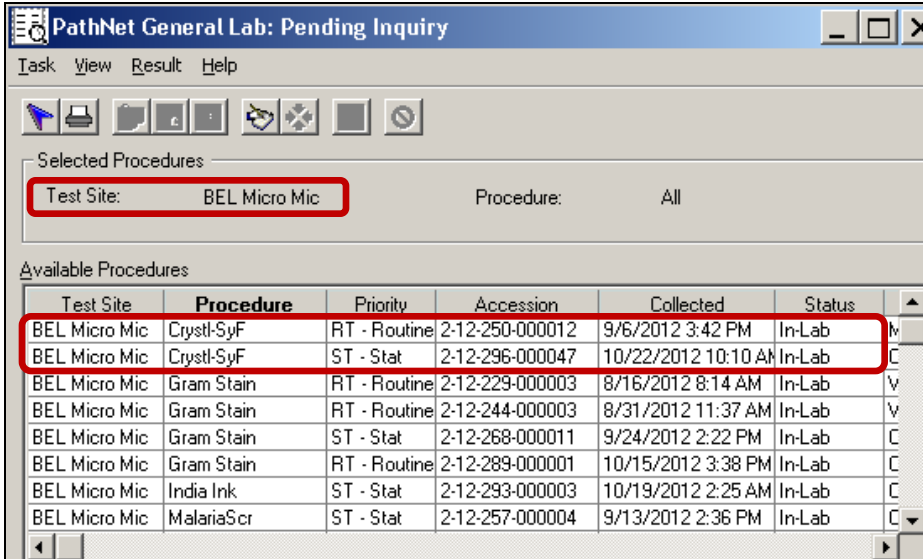
### Changing the Designated Test Site, continued

Step	Action
6	<ul style="list-style-type: none"> <li>Click on <b>T</b>ask on the menu bar.</li> <li>Select <b>Transfer</b> &gt; <b>Selected Orders</b> &gt; <b>All details</b>.</li> </ul> 
7	<ul style="list-style-type: none"> <li>The <b>Transfer</b> window opens.</li> <li>Select the Test Site in the <b>Transfer to Selected Orders To</b> field, clicking on the <b>ellipses</b> as needed.</li> <li>To reprint labels for the specimens being transferred, select <b>Reprint Labels</b>.</li> <li>Click on <b>OK</b>.</li> </ul> 
8	<p>To check if the selected orders are transferred to the new <b>Test Site</b> for resulting:</p> <ul style="list-style-type: none"> <li>Enter new Test Site in the <b>Test Site</b> box, clicking on the <b>ellipses</b> as needed.</li> <li>Click on <b>OK</b>.</li> </ul> 

*Continued on next page*

## Transferring / Tracking Specimens, Continued

### Changing the Designated Test Site, continued


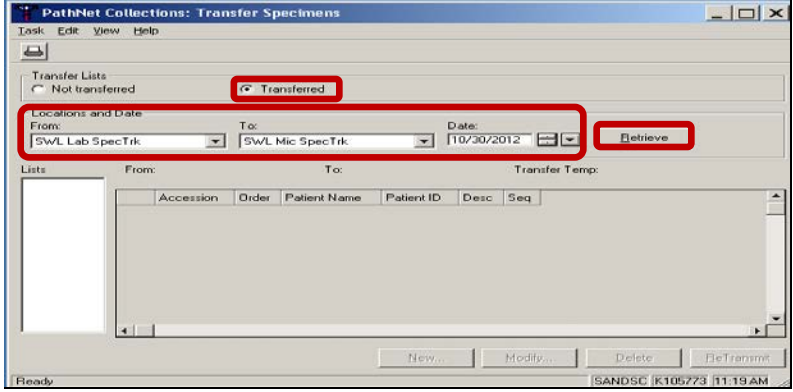
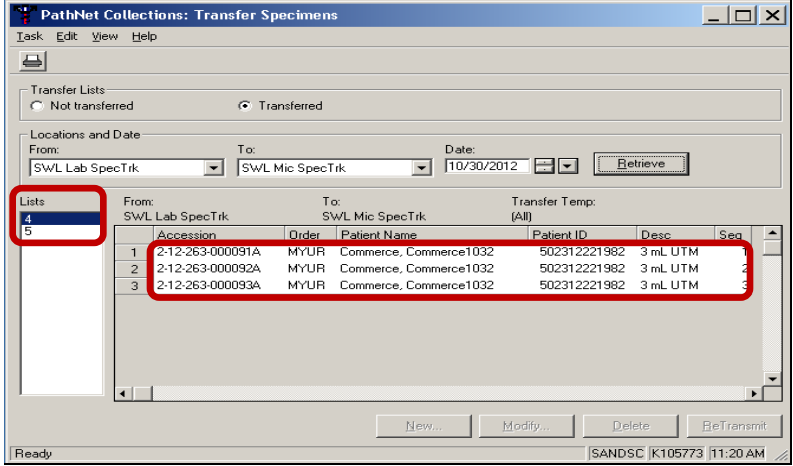
Step	Action																																																						
9	The <b>Pending Inquiry</b> window opens showing the new performing location and the orders that have been transferred.																																																						
 <p>The screenshot shows the 'PathNet General Lab: Pending Inquiry' window. It has a menu bar with 'Task', 'View', 'Result', and 'Help'. Below the menu bar is a toolbar with several icons. The 'Selected Procedures' section shows 'Test Site: BEL Micro Mic' and 'Procedure: All'. The 'Available Procedures' section contains a table with the following data:</p> <table border="1"> <thead> <tr> <th>Test Site</th> <th>Procedure</th> <th>Priority</th> <th>Accession</th> <th>Collected</th> <th>Status</th> </tr> </thead> <tbody> <tr> <td>BEL Micro Mic</td> <td>Crystl-SyF</td> <td>RT - Routine</td> <td>2-12-250-000012</td> <td>9/6/2012 3:42 PM</td> <td>In-Lab</td> </tr> <tr> <td>BEL Micro Mic</td> <td>Crystl-SyF</td> <td>ST - Stat</td> <td>2-12-296-000047</td> <td>10/22/2012 10:10 AM</td> <td>In-Lab</td> </tr> <tr> <td>BEL Micro Mic</td> <td>Gram Stain</td> <td>RT - Routine</td> <td>2-12-229-000003</td> <td>8/16/2012 8:14 AM</td> <td>In-Lab</td> </tr> <tr> <td>BEL Micro Mic</td> <td>Gram Stain</td> <td>RT - Routine</td> <td>2-12-244-000003</td> <td>8/31/2012 11:37 AM</td> <td>In-Lab</td> </tr> <tr> <td>BEL Micro Mic</td> <td>Gram Stain</td> <td>ST - Stat</td> <td>2-12-268-000011</td> <td>9/24/2012 2:22 PM</td> <td>In-Lab</td> </tr> <tr> <td>BEL Micro Mic</td> <td>Gram Stain</td> <td>RT - Routine</td> <td>2-12-289-000001</td> <td>10/15/2012 3:38 PM</td> <td>In-Lab</td> </tr> <tr> <td>BEL Micro Mic</td> <td>India Ink</td> <td>ST - Stat</td> <td>2-12-293-000003</td> <td>10/19/2012 2:25 AM</td> <td>In-Lab</td> </tr> <tr> <td>BEL Micro Mic</td> <td>MalariaScr</td> <td>ST - Stat</td> <td>2-12-257-000004</td> <td>9/13/2012 2:36 PM</td> <td>In-Lab</td> </tr> </tbody> </table>		Test Site	Procedure	Priority	Accession	Collected	Status	BEL Micro Mic	Crystl-SyF	RT - Routine	2-12-250-000012	9/6/2012 3:42 PM	In-Lab	BEL Micro Mic	Crystl-SyF	ST - Stat	2-12-296-000047	10/22/2012 10:10 AM	In-Lab	BEL Micro Mic	Gram Stain	RT - Routine	2-12-229-000003	8/16/2012 8:14 AM	In-Lab	BEL Micro Mic	Gram Stain	RT - Routine	2-12-244-000003	8/31/2012 11:37 AM	In-Lab	BEL Micro Mic	Gram Stain	ST - Stat	2-12-268-000011	9/24/2012 2:22 PM	In-Lab	BEL Micro Mic	Gram Stain	RT - Routine	2-12-289-000001	10/15/2012 3:38 PM	In-Lab	BEL Micro Mic	India Ink	ST - Stat	2-12-293-000003	10/19/2012 2:25 AM	In-Lab	BEL Micro Mic	MalariaScr	ST - Stat	2-12-257-000004	9/13/2012 2:36 PM	In-Lab
Test Site	Procedure	Priority	Accession	Collected	Status																																																		
BEL Micro Mic	Crystl-SyF	RT - Routine	2-12-250-000012	9/6/2012 3:42 PM	In-Lab																																																		
BEL Micro Mic	Crystl-SyF	ST - Stat	2-12-296-000047	10/22/2012 10:10 AM	In-Lab																																																		
BEL Micro Mic	Gram Stain	RT - Routine	2-12-229-000003	8/16/2012 8:14 AM	In-Lab																																																		
BEL Micro Mic	Gram Stain	RT - Routine	2-12-244-000003	8/31/2012 11:37 AM	In-Lab																																																		
BEL Micro Mic	Gram Stain	ST - Stat	2-12-268-000011	9/24/2012 2:22 PM	In-Lab																																																		
BEL Micro Mic	Gram Stain	RT - Routine	2-12-289-000001	10/15/2012 3:38 PM	In-Lab																																																		
BEL Micro Mic	India Ink	ST - Stat	2-12-293-000003	10/19/2012 2:25 AM	In-Lab																																																		
BEL Micro Mic	MalariaScr	ST - Stat	2-12-257-000004	9/13/2012 2:36 PM	In-Lab																																																		

*Continued on next page*

## Transferring / Tracking Specimens, Continued

### Viewing a Transfer List

Follow the steps below to view a transfer list from any current or past date.

Step	Action	Icon																								
1	Click on the <b>Transfer Specimen</b> icon from the Appbar to launch the application.																									
2	<ul style="list-style-type: none"> <li>• The <b>Transfer Specimens</b> window opens.</li> <li>• Select either <b>Transferred</b> or the <b>Not Transferred</b> radio button</li> <li>• Enter the <b>From</b> and <b>To</b> locations, and the <b>Date</b>.</li> <li>• Click on <b>Retrieve</b>.</li> </ul> 																									
3	<ul style="list-style-type: none"> <li>• Multiple transfer list numbers will display in the <b>Lists</b> box.</li> <li>• Highlight the transfer <b>List</b> number to be viewed. The specimens on the selected transfer list displays to the right of the <b>Lists</b> box.</li> </ul>  <table border="1" data-bbox="630 1581 1263 1661"> <thead> <tr> <th>Accession</th> <th>Order</th> <th>Patient Name</th> <th>Patient ID</th> <th>Desc</th> <th>Seq</th> </tr> </thead> <tbody> <tr> <td>2-12-263-000091A</td> <td>MYUR</td> <td>Commerce, Commerce1032</td> <td>502312221982</td> <td>3 mL UTM</td> <td>1</td> </tr> <tr> <td>2-12-263-000092A</td> <td>MYUR</td> <td>Commerce, Commerce1032</td> <td>502312221982</td> <td>3 mL UTM</td> <td>2</td> </tr> <tr> <td>2-12-263-000093A</td> <td>MYUR</td> <td>Commerce, Commerce1032</td> <td>502312221982</td> <td>3 mL UTM</td> <td>3</td> </tr> </tbody> </table>	Accession	Order	Patient Name	Patient ID	Desc	Seq	2-12-263-000091A	MYUR	Commerce, Commerce1032	502312221982	3 mL UTM	1	2-12-263-000092A	MYUR	Commerce, Commerce1032	502312221982	3 mL UTM	2	2-12-263-000093A	MYUR	Commerce, Commerce1032	502312221982	3 mL UTM	3	
Accession	Order	Patient Name	Patient ID	Desc	Seq																					
2-12-263-000091A	MYUR	Commerce, Commerce1032	502312221982	3 mL UTM	1																					
2-12-263-000092A	MYUR	Commerce, Commerce1032	502312221982	3 mL UTM	2																					
2-12-263-000093A	MYUR	Commerce, Commerce1032	502312221982	3 mL UTM	3																					

Continued on next page

## Transferring / Tracking Specimens, Continued

**Non-Controlled Documents** The following non-controlled document supports this procedure.

- Cerner GenLab Participant Guide – Fall 2011 Version 2.1

**Controlled Documents** The following controlled documents support this procedure.

Regional Parent Document Reference Number: SCPMG LIS – 0082 Rev. 2

Job Aid
Create a Transfer List
Modify a Transfer List
Associate a Transfer List to an Outer Tub
Transfer Specimens Using Pending Inquiry
Pocket Guide – Receptionists
Pocket Guide – Lab Assistant
Pocket Guide – Lab Assistant SWL
Pocket Guide – SPD and CSC SWL
Pocket Guide – CLS MC/MOB
Pocket Guide – CLS SWL

<p><b>Author(s)</b> Maureen Ahler, MSQA, MT(ASCP)            Eleonor Arellano, MBA, MBA(HCM), MT(ASCP)            Marian Azuma, CLS            Michelle Borromeo            Karen Brinson, MS, MT(ASCP)            Mario Calantoni, CLS            Cynthia Calderon, SBB, MT(ASCP)            Charles E. Cummings            Ruby Co, CLS            Joshua Evangelista, CLS            Anna Hamilton, MT(ASCP)            Eleanor Ilagan, MT(ASCP)            Julie Kakuda            Allison Kauf</p>	<p>Chongbae Lee, MS, CLS            Betty Lindgren, CLS            Filemon Lizan, CLS            Edna Makalintal, CLS            Linda Mercurio, MAM, MT(ASCP)            Suzanne Nagami            Carmelita Pascua, MT(ASCP)            Chet Phipps, CLS            Stephanie Prien, MS, MT(ASCP)            Carlo Punu, CLS            Alan V. Tran            Diem Trang Vo, CLS            Mary Anne Umekubo, MS, CLS            Janice Wolf, CLS</p>
---	--