



KAISER PERMANENTE®

DOCUMENT NUMBER:
DOCUMENT TITLE:
DOCUMENT NOTES:

LOCATION:	VERSION:
DOC TYPE:	STATUS:

EFFECTIVE DATE:	NEXT REVIEW DATE:
RELEASE DATE:	EXPIRATION DATE:

AUTHOR:	PREVIOUS NUMBER:
OWNER:	CHANGE NUMBER:

Suspension of Autoverification for IRIS IQ200 and Remisol

**Purpose or
Principle or
Introduction**

This procedure provides instructions on how to suspend and re-enable autovalidation or autoverification in Remisol Advance for Chemistry, ImmunoChemistry, Coagulation and Iris IQ 200 Urinalysis System.

Scope

All CLS and MLT who work on systems with autoverification capabilities have the access and right for this document usage.

Policy

Suspend autoverification if:

- a. QC results are unacceptable even after thorough investigation and corrective action.
- b. Instrument malfunction.
- c. Failed calibration.
- d. Reagent problem.
- e. Any other situation where instrument results are suspected to be adversely affected, e.g. high temperature in the lab.

Re-enable autoverification if:

- a) QC results are acceptable.
- b) Instrument problem has been corrected.
- c) Calibration has passed.
- d) Reagent problem has been solved.
- e) Any other situation has been corrected and instrument result are now acceptable.

Note:

If unsure of what to do, it is good laboratory practice to find policies/instructions on Master Control, consult Manager, lead CLS or call the vendor hotline.

Continued on next page

Suspension of Autoverification for IRIS IQ200 and Remisol, Continued

IRIS IQ200

**Suspension of
 autoverification**

Step	Action
1.	To suspend autovalidation on IQ200 <ul style="list-style-type: none"> ▪ Log on to IQ200 computer as Manager ▪ Click on Instrument ▪ Go off line Click on Settings ▪ Click on Auto-Release ▪ Uncheck Enable Auto- Release ▪ Click on OK
2.	Verify that autoverification is successfully turned off by running a previously known negative patient accessioned with a non-patient Health Connect test MRN and verify that results are not auto released. Document on daily checklist.

IRIS IQ200

**Re-enable of
 autoverification**


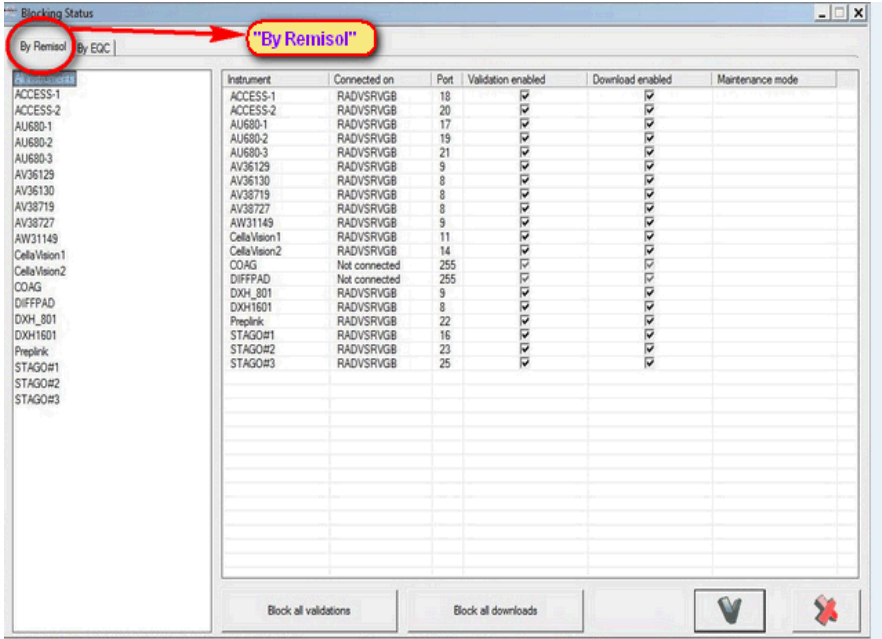
Step	Action
1.	To re-initiate autoverification on IQ200 <ul style="list-style-type: none"> ▪ Log on to IQ200 computer as manager ▪ Click on Instrument ▪ Go off line ▪ Click on Settings ▪ Click on Auto-Release ▪ Check the enable Auto-Release box ▪ Click on OK
2.	Verify that autoverification is successfully turned on by running a previously known negative patient accessioned with a non-patient Health Connect test MRN and verify that results are not auto released. Document on daily checklist.

Continued on next page

Suspension of Autoverification for IRIS IQ200 and Remisol, Continued

REMISOL

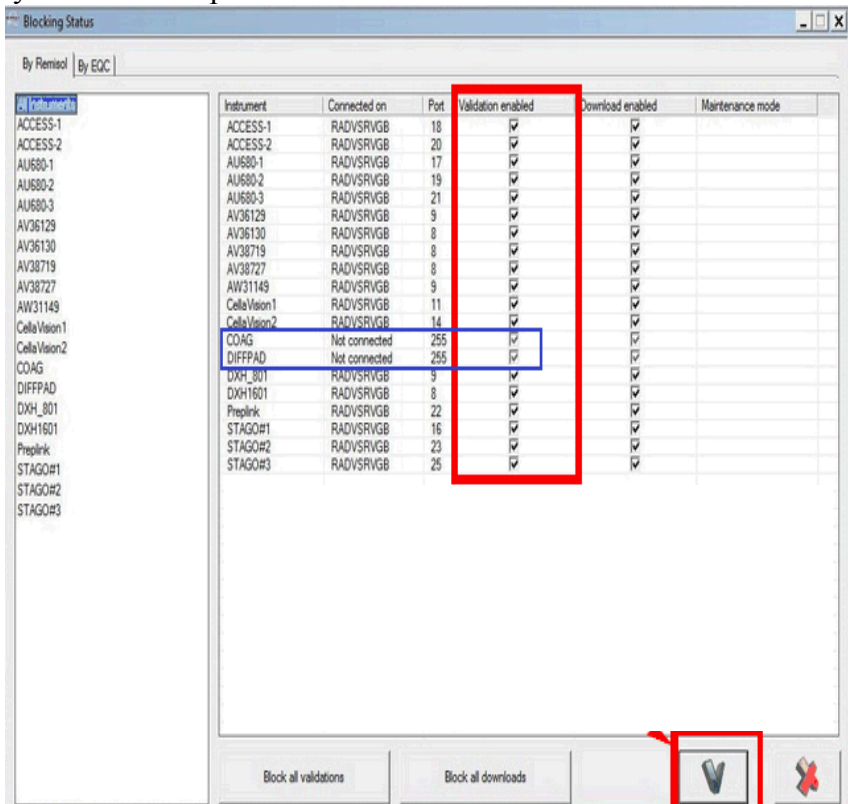

**Suspension of
 autoverification**

Steps:	Actions:
Note	Use the Manager’s Remisol station in room # 1835.
1	Click the STOP icon located in the Tool Bar, 
2	Select the “By Remisol” tab located in the upper left corner of the <i>Blocking Status</i> box. 


Continued on next page

Suspension of Autoverification for IRIS IQ200 and Remisol, Continued

Continuation
REMISOL
 Suspension of
 autoverification

Steps:	Actions:
3	<p>In the “Validation Enabled” column, unmark the box of the instrument you wish to suspend autoverification.</p> 
4	Save by clicking the Green V button.
5	<p>This RED HAND will now display on the “STOP” icon, .</p>

Suspension of Autoverification for IRIS IQ200 and Remisol, Continued

	Steps	Actions
REMISOL Re-enable autoverification	Note	Re-enable of autoverification is only for Managers or Lead CLS. Use the Remisol in the Manager’s office.
	1	Log-in with Administrator’s username and password.
	2.	Click the Stop icon, 
	3.	Select the “By Remisol” tab located on the upper left corner of the Blocking Status box.
	4.	Check the boxes under the Validation Enabled column to re-enable autoverification.
	5.	Click the green V to save changes. The Stop icon without the red hand will now display on the Remisol’s Toolbar.
	6.	Log-out of the Remisol by pressing together CTRL + B.
	7.	Verify that all Remisol stations took in the re-enabled autovalidation.

**Safety or
 Special Safety
 Precautions**

Refer to the safety manual for general safety requirements.

Definitions

Autoverification or Autovalidation	Results which are within the normal ranges are auto validated or autoverified if it meets pre-defined criteria.
------------------------------------------	-----------------------------------------------------------------------------------------------------------------

Author(s) Anita Jo-Ann Calalang, CLS

Signature Manifest

Document Number: LAMC-PPP-0775

Revision: 01

Title: Suspension of Autoverification for IRIS IQ200 and Remisol

All dates and times are in Pacific Standard Time.

Suspension of Autoverification

Change Request

Name/Signature	Title	Date	Meaning/Reason
Anita Calalang (O880586)	LAB INFO SYS APPLIC SPECIALIST	04 Aug 2017, 02:32:59 PM	Approved

Quality Approval

Name/Signature	Title	Date	Meaning/Reason
Jocelyn Javier (T684676)	Assist. ADA	09 Aug 2017, 11:32:17 AM	Approved

Operations Approval

Name/Signature	Title	Date	Meaning/Reason
Julie Toti (K084521)	DIR AREA LAB	10 Aug 2017, 10:47:02 AM	Approved

Final Approval

Name/Signature	Title	Date	Meaning/Reason
Hedyeh Shafi (I086749)	Pathologist	14 Aug 2017, 11:34:15 AM	Approved

Set Effective Date

Name/Signature	Title	Date	Meaning/Reason
Anita Calalang (O880586)	LAB INFO SYS APPLIC SPECIALIST	25 Sep 2017, 01:13:37 PM	Approved

Quick Approval

Approve Now

Name/Signature	Title	Date	Meaning/Reason
Anita Calalang (O880586)	LAB INFO SYS APPLIC SPECIALIST	03 Oct 2017, 11:47:39 AM	Approved