


Table of Contents	Day Of Collection Workflow	Introductory Information	Procedures	Packaging	Resources	Version History
			Clinical Trial	NanoCool		
			Commercial	Credo Cube		
				QRG		
				Training		
				Deviations		
				Definitions		
				Attachments		

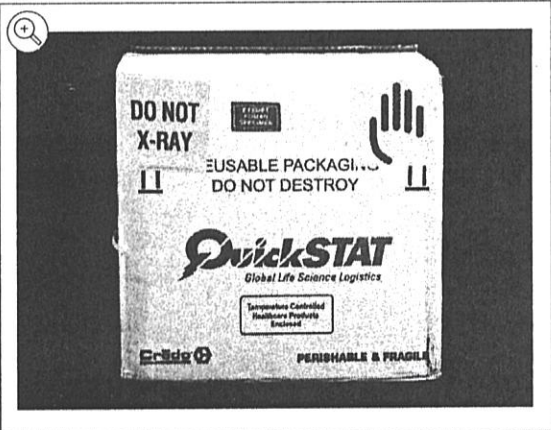


Label the outer cardboard corrugate the Credo arrived in with the following labels:

- ‘DO NOT X-RAY’
- ‘Exempt Human Specimen’
- BMS logo


Step 4

COI Verification: Prior to affixing the Shipping Address Label, verify the JOIN on the MNC Bag Label, Shipping Address Label, and CSMCC* matches for the product being packaged. Affix the verified Shipping Address label to the side of the outer cardboard box.

** Or the MNC Collection Procedure Record if using the manual backup procedure and the CSMCC is not available.*

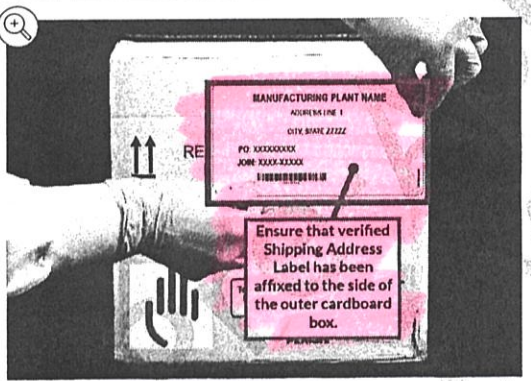

Ensure that the verified Shipping Address Label has been affixed to the side of the outer cardboard box.

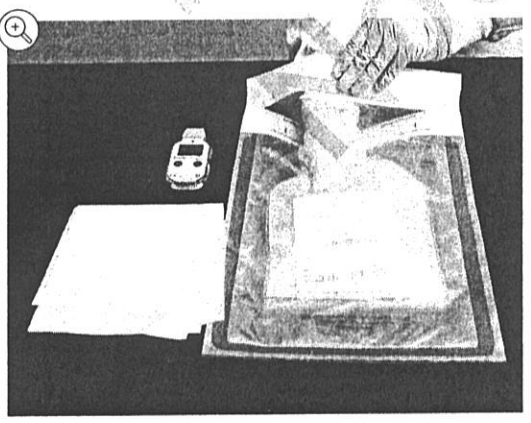

Place the MNC Product and absorbent sheet into the specimen transport bag with the MNC Bag label facing outward.

Note: MNC Bag label should be visible and not obstructed. The label should be facing the side of specimen transport bag without the biohazard print, so the label is visible while inside the bag.

Expel the air from the specimen transport bag, remove the adhesive, and seal the bag.