

Change Summary: Updated Shipper for BMS Cell Therapy Drug Products

BMS Confidential Information. © 2026 Bristol-Myers Squibb Company.
03/2026 HE-US-2600245

Disclaimers

- This presentation is intended for training on operational matters for implementing CAR T Cell therapy at Sites of Care.
- Please note the intent is to discuss operational considerations and not to promote any products.
- The content of this presentation and related conversations are subject to a confidentiality agreement and should not be disseminated to anyone not covered by the confidentiality agreement.

Change Summary

- BMS is implementing the Cryoport HV3 General Purpose Shipping System, featuring a new exterior shipping case for cryogenic shippers
- There is no change to:
 - The interior shipper chamber which holds drug product
 - Thermal performance or validation of the cryogenic shipper
 - Data logger or temperature monitoring process
 - Product handling procedures
- Change will be implemented starting 1 Apr 2026

Current shipper:

- Cylindrical outer shell with rounded top
- Shipping container requires separate dolly or mechanism for transport
- Hanging document pouches inside
- Paper instructions included
- Cannot be stacked or palletized



Updated Shipper (as of April 1, 2026)

- Cube-shaped outer shell with built-in dolly system
- Integrated document compartment (no hanging pouches)
- Designated exterior label panel
- Stackable design for transport
- Same interior chamber with 8-day validated hold time at cryogenic temperatures as current shipper

Exterior Shipper Specifications

Product Specifications	Updated shipper as of April 1, 2026- HV3 Shipping System	Current shipper- HV2 Shipping System
Image		
Temperature Range	-150C or colder	-150C or colder
Dynamic Hold Time	Up to 8 days	Up to 8 days
Exterior Dimensions	22" L x 22" W x 28" H 558.8 mm x 558.8 mm x 711.2 mm	22" D x 28" H 558.8 mm x 711.2 mm
Interior Chamber Dimensions	8.5" (215.9 mm) diameter 12" (304.8 mm) height Max commodity height 11.5"	8.5" (215.9 mm) diameter 12" (304.8 mm) height Max commodity height 11.5"
Dry Weight	Up to 73 lb / 33 kg	Up to 57lb / 26 kg
Charged Weight	Up to 92 lb / 42 kg	Up to 77 lb / 35 kg


Exterior Images



Exterior Images

Waybills & shipping labels

Dolly handle

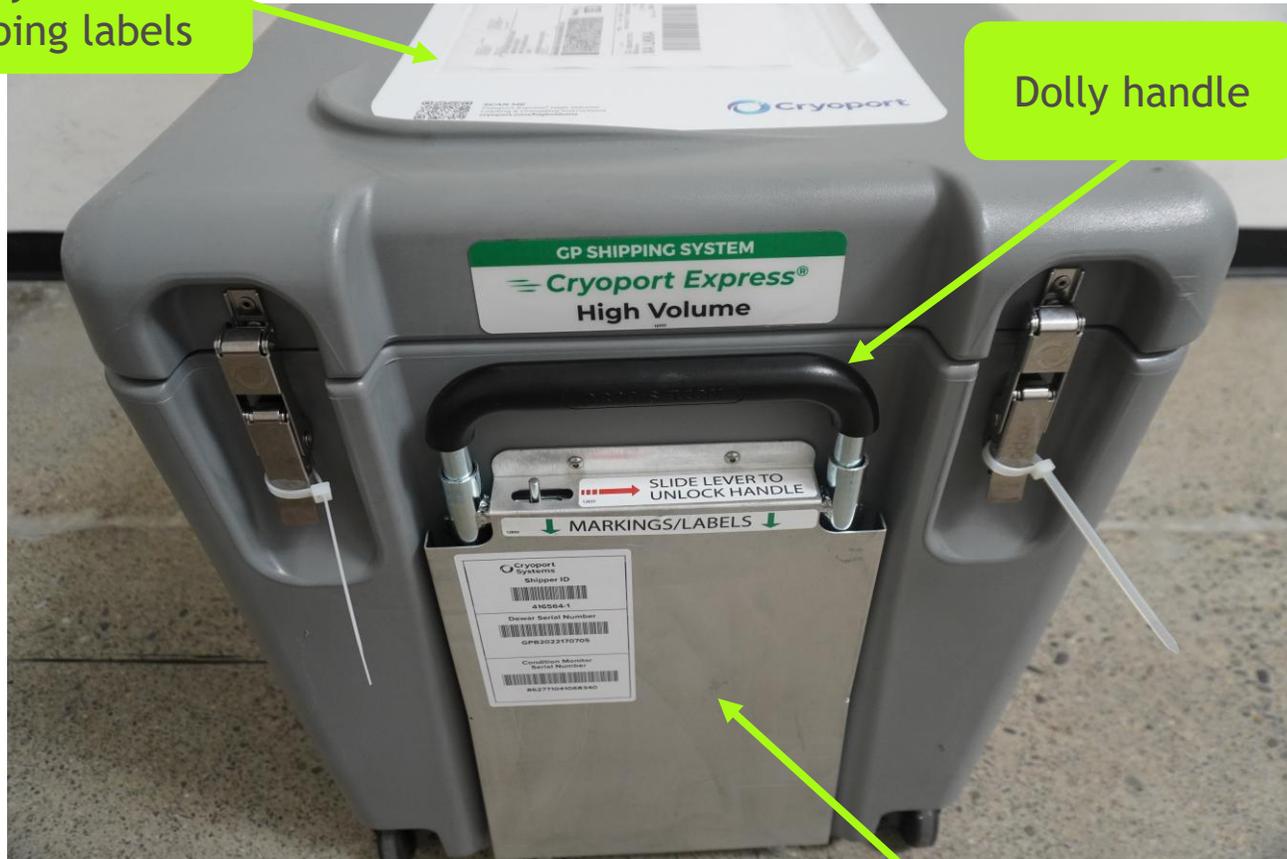

Shipper expiry label panel



Zip tie and latch closure

Opening the shipper

1 Cut the zip-tie on each latch



2 Depress the button on the underside of the latch and lift to release




Push up button
on the underside
of the latch



Document layout change - inside

Current

There Will Be Two Shipping Pouches Inside the Cryoshipper

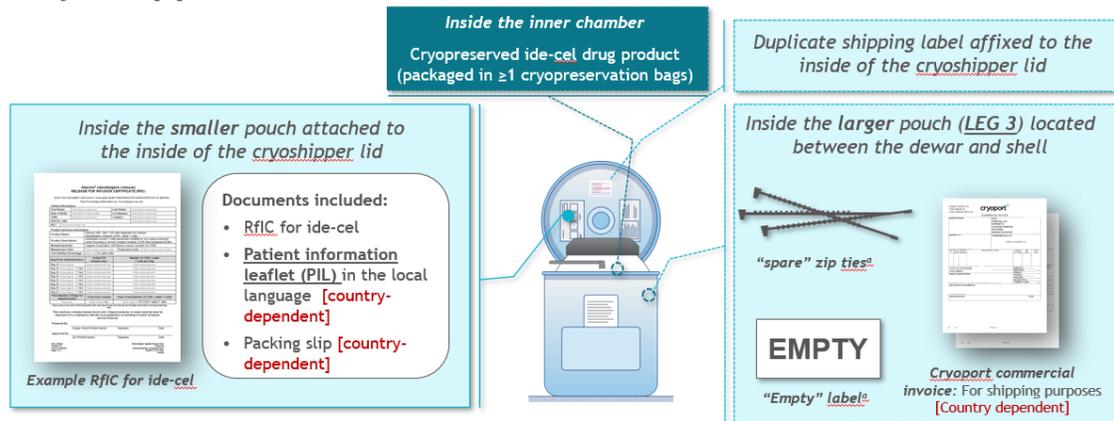

¹Treatment sites may decide whether to use the label, or zip ties. If the site does not need to use these items, the site may leave them untouched in the LEG 3 pouch.

PIL, patient information leaflet; RfIC, Release for Infusion Certificate.
References: 1. BMS internal files. 2022. 2. Abocma, idocabazone vicleucel) Release for Infusion Certificate. Varies from market to market.

Contact BMS at [PHONE NUMBER] with questions.

Bristol Myers Squibb | ide-cel Clinician Handling Guide

Subject to confidentiality agreement. Do not disseminate without prior authorization by BMS. | 28


Future



2 plastic pouches included in the document compartment:

- Leg 2 will contain:
 - Release for Infusion Certificate
 - Patient information leaflet [country dependent]
 - Packing slip [country dependent]
 - Secondary BMS shipper expiry label**
- Leg 3 will contain:
 - Empty label
 - Spare zip ties
 - Commercial invoice [country dependent]

Accessing the Document Compartment

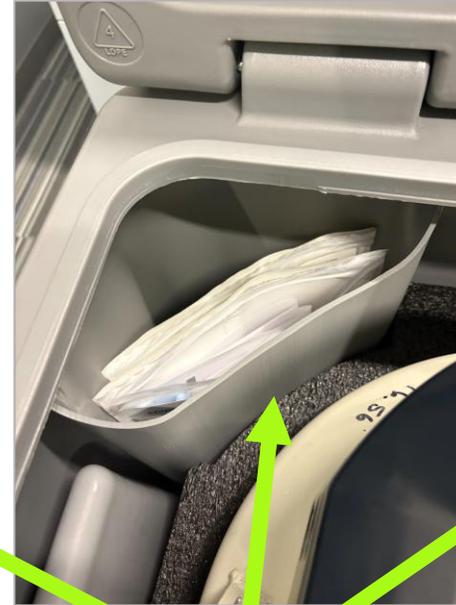

WARNING: The lid opens and rests at a 90-degree angle. Close it carefully to avoid slamming, as pushing it will cause it to shut abruptly.

Document compartment is clearly visible and accessible when the exterior lid is open

Using the Built in Dolly to Move the Shipper

1

Slide lever to unlock handle and pull handle up with your other hand.



2

Lift handle until you hear an audible click. Handle can extend to two different heights. Tilt the shipper and use wheels to pull like a suitcase.



Label Placement

Label requirements may vary by country





UN3245 sticker

Drug name: Bristol Myers Squibb
 Transporter used No.
 Full Delivery address for education:
 -Site name
 -Street
 -Town
 -Postcode
 Packaging Expiry date/time:

Shipper expiry label



Do Not X-Ray sticker
[country dependent]

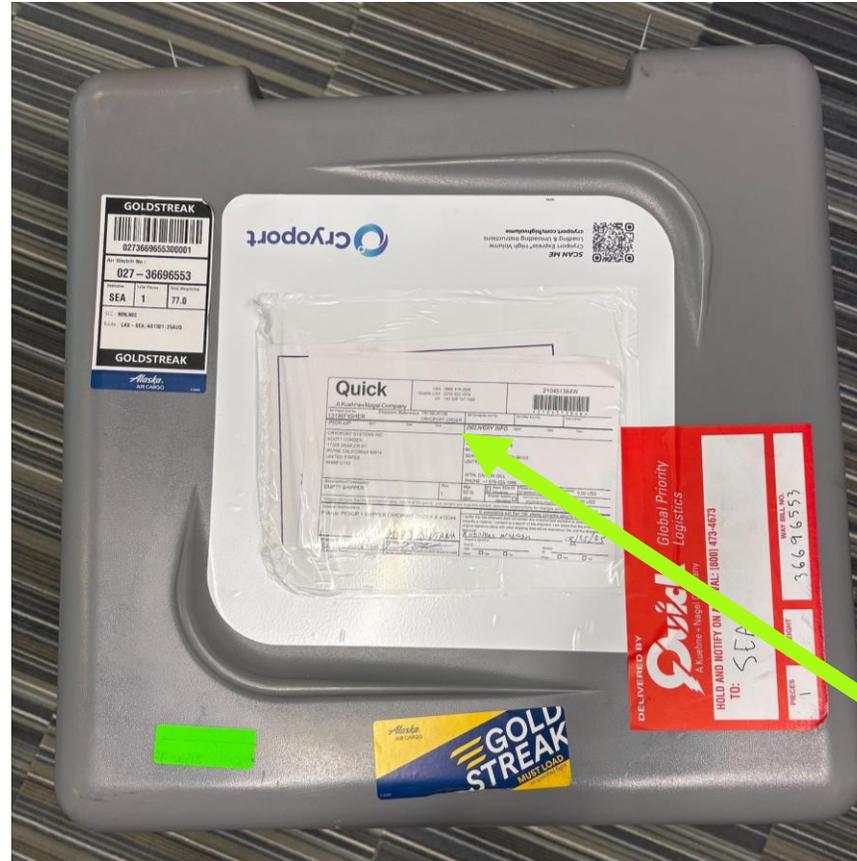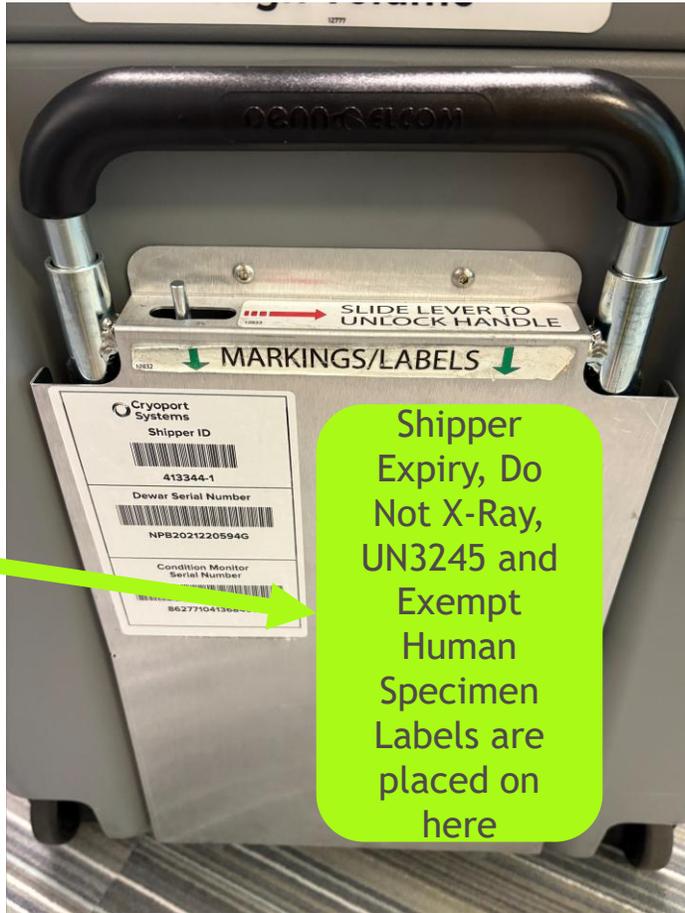

MARKEN
PACKAGE # 1 of 2

64610830525

PLACEMENT

Package to Place
Type: No. of Packages: No. of Pieces:
 Shipper: Receiver: Date:

Event Job Number: 04610830525

Special Reporting Instructions

Consignee Signature
Name: Title:
 Received in Good Order and Condition

Waybill

Waybill and return label [country dependent] are placed here

Return Shipper Instructions: Secure the latch with zipties



Thread the zip tie through the hole and close it around the latch




Follow BMS Handling Instructions



ALWAYS FOLLOW
BMS INSTRUCTIONS
FOR CELL THERAPY
DRUG PRODUCT AND
SHIPPER HANDLING.

DISREGARD
CRYOPORT.COM
INSTRUCTIONS
LINKED TO AT THIS
QR CODE.

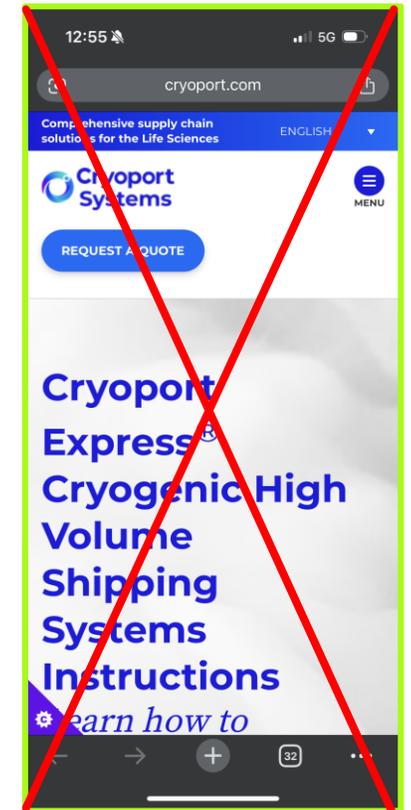

Contact BMS CT360 Scheduling for any questions

Scheduling@celltherapy360.com