



SCC-LIS_TIP OF THE WEEK

MODULE: SoftLab/SoftMic	SOFTWARE VERSION: 4.0.8
TOPIC: Instrument Menu Test Rerun	ISSUE DATE: 19APRIL2017
DISTRIBUTION: End Users_All	PAGE: 1 of 1
ISSUED BY: Joel Droogers Territorial LIS Administrator	

TOPIC/QUESTION:

How do I get my analyzer to rerun tests without having to manually program?

ANSWER/TIP:

Open Instrument Menu:

- Rerun one test/selected tests:
 - Click on the test to rerun – then click Rerun→this will send the test back to the loadlist
 - Repeat for each test you wish to rerun

Patient Info
 Name (last, first, middle): TESTING, STANTON
 CLIENT S:00000685 Sex: F Age: 116 Ward: YEL Room:
 Order Info
 Order: 72010162-04 Prior: R Seq: 0075 Spec: SST Collected: 01/12/2016 11:00
 Tests: CA, K

Itst	T Name	Result	R	Flags	M	Prev Res	Prev Date	Error	Status	Tech	DateV
1	308 Potassium	4.6				4.5	01/12/2016		C		
2	318 Calcium	2.16				2.50	14/07/2016		C		
3	950 Hemolysis Check, Vitros	<15							5 C		
4	952 Turbidity Check, Vitros	<20							5 C		
5	951 Icterus Check, Vitros	<2							5 C		

- Rerun EVERY test on the order:
 - Click on the first test in the list
 - Hold the shift key and use the down arrow to highlight all the tests
 - Hit the Rerun button→this will send all the tests highlighted back to the loadlist

Patient Info
 Name (last, first, middle): TESTING, STANTON
 CLIENT S:00000685 Sex: F Age: 116 Ward: YEL Room:
 Order Info
 Order: 72010162-04 Prior: R Seq: 0075 Spec: SST Collected: 01/12/2016 11:00
 Tests: CA, K

Itst	T Name	Result	R	Flags	M	Prev Res	Prev Date	Error	Status	Tech	DateV
1	308 Potassium	4.6				4.5	01/12/2016		C		
2	318 Calcium	2.16				2.50	14/07/2016		C		
3	950 Hemolysis Check, Vitros	<15							5 C		
4	952 Turbidity Check, Vitros	<20							5 C		
5	951 Icterus Check, Vitros	<2							5 C		

- Load the sample on the analyzer→rerun results will append in Instrument Menu

72010162-04	C	0075	5	11:30	CK, CCA, CHEM, CCT, CTUR
72010162-04	P	0077	5	11:39	PK, PCA, CHEM, CCT, CTUR