



SCC-LIS_TIP OF THE WEEK

MODULE: SoftLab/SoftBank	SOFTWARE VERSION: 4.0.8
TOPIC: Blood Bank History Check	ISSUE DATE: 02FEB2018
DISTRIBUTION: End Users_All	PAGE: 1 of 1
ISSUED BY: April Darrach, Territorial LIS Administrator	

TOPIC/QUESTION:


What is this HXCK test that keeps reflexing whenever I order a Type and Screen (TS)?

ANSWER/TIP:

HXCK = Blood Bank History Check

To ensure that we are recording relevant blood bank patient history – we have created a HXCK test code that includes: checking the card file, checking for previous transfusions and pregnancies. These tests are resulted in the SoftBank module. This test code, once LIVE, will be specific to Fort Smith, Inuvik and Stanton sites.

EXAMPLE:

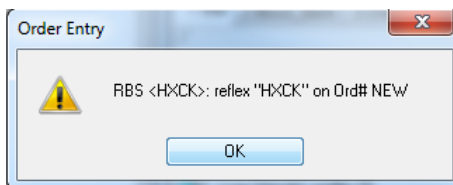

SoftLab – results are auto-verified with the result of “See Blood Bank”

T Ord	T Name	Result	FI	Flags	M	Prev Res	Prev Date	C	Status	Tech	DateV	TimeV
1	HXCK Card File Read	See Blood Bank							LC V		2018-01-27	13:53
2	HXCK Transfusion within the last 3 months?	See Blood Bank							LC V		2018-01-27	13:53
3	HXCK Pregnancy within the last 3 months?	See Blood Bank							LC V		2018-01-27	13:53

SoftBank – results must be entered for each individual test. They are resulted with “Yes”, “No” or “UA”(unavailable). See Transfusion Medicine Manual for further information

History Check		
Card File Read?		requested
Previous Tx?		requested
Prev. Pregnancy?		requested