

# NWT LIS TIP OF THE WEEK

**Module:** SoftLabBank

**Software Version:** 4.0.8

**Topic:** Closing BBID Properly

**Issued On:** 22NOV2018

**Distribution:** All Users

**Page:** 1 of 1

**Issued By:** M-L Dufresne, Territorial LIS Administrators

## TOPIC/QUESTION:

What should I do when the BBID window pops up for tests where a blood bank arm band is not applicable?


## ANSWER/TIP :

Leave the BBID field blank and hit "OK".

Avoid cancelling or "X-ing" out of the window. If you do this, the collection information will not save.


## EXAMPLE :

### Correct

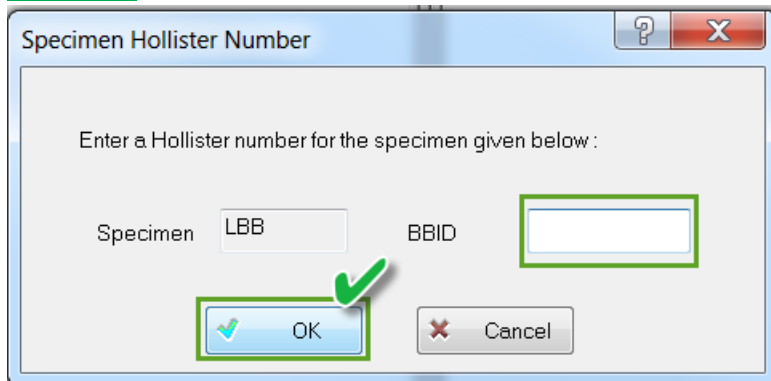

The screenshot shows a dialog box titled "Specimen Hollister Number". It contains the text "Enter a Hollister number for the specimen given below :". Below this, there are two input fields: "Specimen" with the value "LBB" and "BBID" which is empty. The "BBID" field is highlighted with a green border. At the bottom, there are two buttons: "OK" and "Cancel". The "OK" button is highlighted with a green border and has a green checkmark above it, indicating the correct action.

### Incorrect

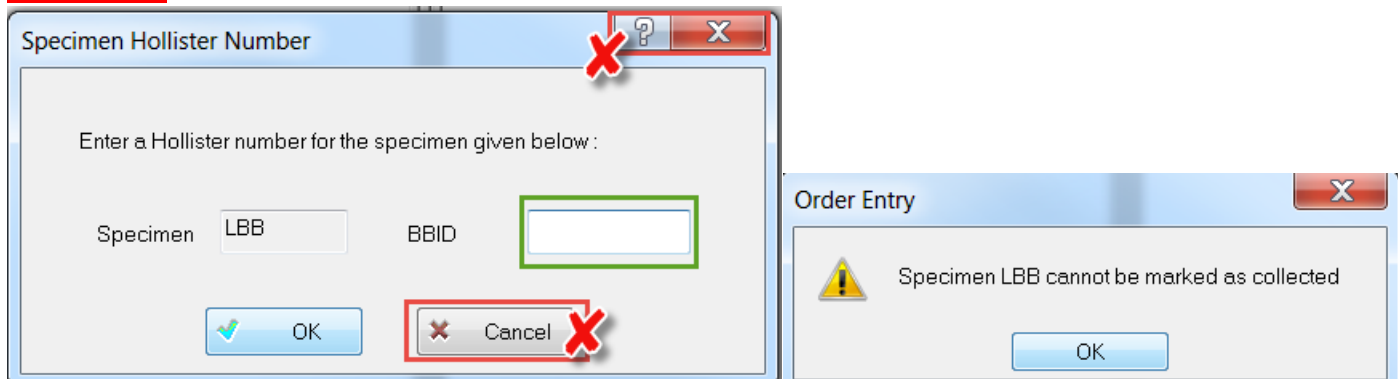

The screenshot shows the same "Specimen Hollister Number" dialog box as above, but with a red 'X' over the "Cancel" button, indicating that clicking it is incorrect. To the right, a second dialog box titled "Order Entry" is shown, displaying a warning icon and the message "Specimen LBB cannot be marked as collected". This dialog box has an "OK" button at the bottom.