

NWT LIS TIP OF THE WEEK

Module: SoftLab	Software Version: 4.0.8
Topic: Renal Panels_SLAB	Issued On:11JUNE2019
Distribution: All Users_SLAB	Page: 1 of 1
Issued By: M-L Dufresne, Territorial LIS Administrator	


TOPIC/QUESTION:

I'm trying to accession an order on a renal patient that has upwards of 50 tests on it!!!! How can I best manage this?

ANSWER/TIP:

LIS orders can have a **maximum** of **60** tests on them. It is ill-advised to accession this many tests on an order as this will potentially not leave any room on the order for reflexed testing that may occur once samples are run.

A special ordering keypad has been created for renal panels that should facilitate ordering the in-house tests on one order number and the referral tests on a separate order number. This will avoid having one order number with too many tests on it.

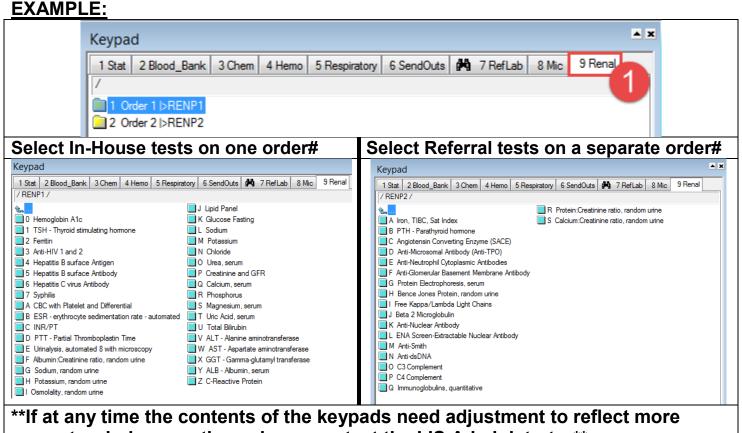

current ordering practises please contact the LIS Administrator**