

NWT LIS TIP OF THE WEEK

Module: SoftLab/Mic

Software Version: 4.0.8

Topic: RSV Testing in Core Lab

Issued On: 17Sep2019

Distribution: Stanton Users

Page: 1 of 1

Issued By: M-L Dufresne, Territorial LIS Administrator

TOPIC/QUESTION:

Now that core lab is doing RSV testing, what do I need to do in the LIS?

ANSWER/TIP:

Test	RSV
Resulting Worklist Template	SCRN
TQC Lot Record ID	RSV
TQC Frequency	Per Box (Create Ad-Hoc Order)
RBS Rule to add Viral Culture (VIRC)	@VIRC
Scheduled Go-Live Date	12Sept2019
Sharepoint Task	01396

EXAMPLE:

Order Test Code "RSV"

"VIRC" will reflex by RBS Rule

Go to SCRN resulting Worklist

Use keypad to record observations from RSV testing card. Based on observations, system will interpret result and add appropriate result comment. Only the result comment will appear on the report (observations will not)

@RSVI	@RSVP	@RSVN
INVALID testing results obtained by immunochromatographic assay Sent to Provincial Laboratory for Public Health for viral culture	POSITIVE for RSV antigen by immunochromatographic assay Sent to Provincial Laboratory for Public Health for confirmation by viral culture	NEGATIVE for RSV antigen by immunochromatographic assay Sent to Provincial Laboratory for Public Health for confirmation by viral culture