



Stanton Territorial Hospital
P.O. Box 10, 550 Byrne Road
YELLOWKNIFE NT X1A 2N1

Document Name:
Dispensette Sterility Testing Job Aid

Document Number:
MIC70210

Distribution:
Microbiology Instrumentation Manual

Date Issued:
19 March 2021

Order standing order in LIS:

- Patient Last Name: **TEST**
- Patient First Name: **MICRO STERILITY**

- **CXENV** will be ordered
- **Collect, receive and plate**
- **MICRO OE Comments:** Vitek Dispensette

Label 2 THIO broths with specimen labels

Dispense 3mL THIO broth

Read broth daily for 3 days

THIO broth is cloudy (indicates growth)

Record: **"Cloudy-AUTOCLAVE dispenser ^CLAVE"**

Report: **"Cloudy-AUTOCLAVE dispenser"** and finalize

Discard saline bottle and autoclave dispensette.

THIO broth is clear (indicates no growth)

Record: **"No Growth"** in LIS daily for 3 days

Report: **"No Growth After 3 Days"** and finalize