


NWT LIS TIP OF THE WEEK

Module: SoftLabMic	Software Version: 4.0.8
Topic: Barcode Scanners	Issued On: 2023-11-24
Distribution: End All_Users	Page: 1 of 1
Issued By: Lena Schofield, Territorial LIS Administrator	

TOPIC/QUESTION: We received a new barcode scanner but it doesn't scan specimens the same as the old one.

ANSWER/TIP: The Honeywell 1470g Barcode Scanners have slightly different settings than the old scanners, here are the new barcodes to scan to reset then apply new settings. Scan one by one and wait for the beeps that the confirm the setting (some sound different).

The following barcode defaults all of the work groups to the factory default settings.

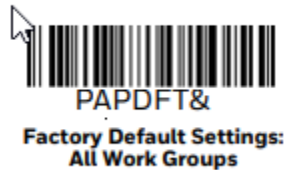

Scanning this barcode reboots the scanner and causes it to relink with the base or Access Point.

