

TRAINING UPDATE

Lab Location: WAH
Department: Micro

Date Distributed: 11/4/2013
Due Date: 11/30/2013
Implementation: 12/1/2013

DESCRIPTION OF PROCEDURE REVISION

Name of procedure:
Microbiology 35C CO2 Incubator log, AG.F19.3
Description of change(s):
Revise to be WAH site-specific log Change to record readings to daily (instead of each shift) This revised log will be implemented on December 1, 2013

Document your compliance with this training update by taking the quiz in the MTS system.



MONTH / YEAR _____

Temperature, humidity and carbon dioxide are documented each day. The carbon dioxide percentage is checked with the Fyrite monthly. The CO₂ gauge is monitored daily. Change tank when gauge indicates <100 PSI. **Never** wait until the CO₂ tank is empty.

When the temperature is higher or lower than the acceptable range, turn the thermostatic control up or down slightly as needed. Acceptable temperature range is 33-37°C.

Any adjustments made should be documented on this sheet and the lead tech or supervisor notified.

Date	Fyrite 4.5-7.5% (record monthly)	Comments	Tech

Date	Temp 33-37°C	Humidity Water in pan (Y/N)	%CO ₂ 5-7%	CO ₂ pressure change tank at <100	Tech
1					
2					
3					
4					
5					
6					
7					
8					
9					
10					
11					
12					
13					
14					
15					
16					
17					
18					
19					
20					
21					
22					
23					
24					
25					
26					
27					
28					
29					
30					
31					

Weekly review:	Weekly review:	Weekly review:
Weekly review:	Weekly review:	Monthly review: