

TRAINING UPDATE

Lab Location: SGMC & WAH
Department: Core

Date Distributed: 7/29/2015
Due Date: 8/19/2015
Implementation: 8/19/2015


DESCRIPTION

Name of procedure:
How to Result a Urine Culture Job Aid AG.F330.0
Description:
<p>This is a new “form” created to assist staff on the proper keys to use when report urine cultures.</p> <p>This Document will be implemented on August 19, 2015</p>

Document your compliance with this training update by taking the quiz in the MTS system.


Sterile sources/specimen description (SDES) - 4 plates and held for 2 days	
UBLAD	BLADDER URINE
UINOUC	URINE IN/OUT CATH
UKID	URINE, KIDNEY
URCYS	URINE CYSTOSCOPY
USUPR	URINE, SUPRAPUBIC

Non sterile sources/specimen description (SDES) - 2 plates and held for 1 day	
Any urine source NOT mentioned above, <i>Example</i> URCC – Urine clean catch	

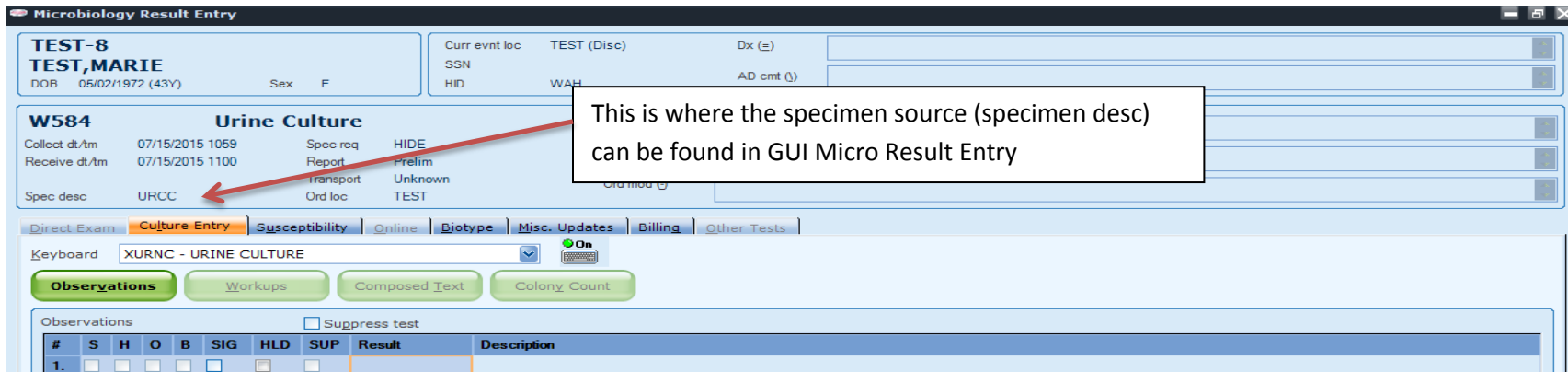

This is where the specimen source (specimen desc) can be found in GUI Micro Result Entry

Resulting (F8 displays keyboard in Micro Result Entry; F8 again will turn if off)

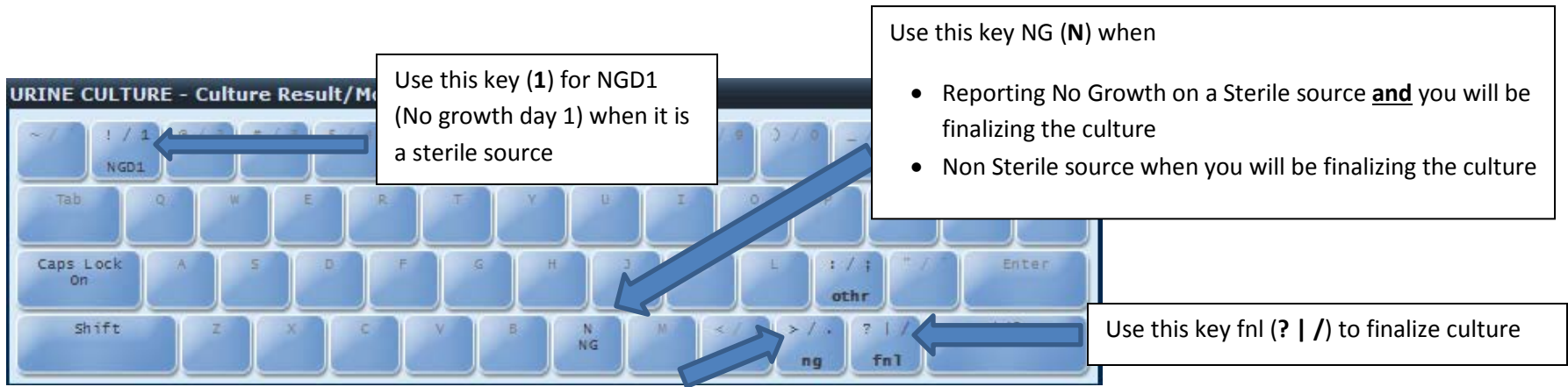

DO NOT USE THIS KEY ng (> / .) TO RESULT