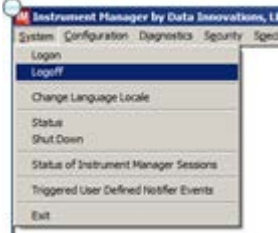
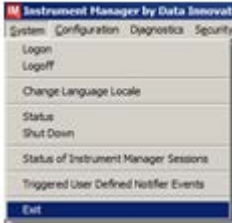



Date: May 30, 2017

INTERNAL LABORATORY ALERT

Subject: Logging out of DI (Instrument Manager)

The Laboratory is informing you of the following:

<p>Logging out of DI (Instrument Manager)</p>	<p>Reminder that logging out of DI is two-step process. If you only perform 1 of the 2 steps, your computer is still tied to DI.</p>
<p>Step 1: Log off</p>	<p>Select System from the upper left corner - Point and Click Pop-up message displays confirming that you want to logoff. Click on Yes.</p> 
<p>Step 2: Close DI application</p>	<p>Two options for closing DI</p> <ol style="list-style-type: none"> Select Exit from the upper left corner - Point and Click. DI application will close.  <ol style="list-style-type: none"> From the main screen, click on the X in the upper left corner. 
<p>Msabonis/053017</p>	