

FREEZER PM LOG – JEWETT JPL1230A

St. Joseph Medical Center Tacoma, WA
 St. Clare Hospital Lakewood, WA
 St. Elizabeth Hospital Enumclaw, WA
 St. Francis Hospital Federal Way, WA
 St. Anthony Hospital Gig Harbor, WA
 Highline Medical Center Burien, WA
 PSC

Year _____ Site: **SAH** Jewett Freezer **JPL1230A** BioMed # **A12568**

Temperature alarm must activate before rising to -18 C.
Backup batteries must be replaced after 1 year.

QUARTERLY PM:

Task	1 st Quarter		2 nd Quarter		3 rd Quarter		4 th Quarter	
Clean	Performed?		Performed?		Performed?		Performed?	
Condenser grill	Y N		Y N		Y N		Y N	
Check	OK?	Replaced ?	OK?	Replaced ?	OK?	Replaced ?	OK?	Replaced ?
Backup battery for built-in chart recorder	Y N	Y N	Y N	Y N	Y N	Y N	Y N	Y N
Backup battery for monitoring system	Y N	Y N	Y N	Y N	Y N	Y N	Y N	Y N
Test	Alarms?	Clears?	Alarms?	Clears?	Alarms?	Clears?	Alarms?	Clears?
Power failure alarm	Y N	Able to clear manually? Y N	Y N	Able to clear manually? Y N	Y N	Able to clear manually? Y N	Y N	Able to clear manually? Y N
Door open alarm	Y N	Y N	Y N	Y N	Y N	Y N	Y N	Y N
Tech ID								
Date								
Reviewed by								
Date								

Task	1 st Quarter		2 nd Quarter		3 rd Quarter		4 th Quarter	
Check probe temperature calibration for monitoring system	Temp Therm _____C Temp digital _____C OK? Y N Readout recalibrated? Y NA		Temp Therm _____C Temp digital _____C OK? Y N Readout recalibrated? Y NA		Temp Therm _____C Temp digital _____C OK? Y N Readout recalibrated? Y NA		Temp Therm _____C Temp digital _____C OK? Y N Readout recalibrated? Y NA	
Alarm Check <i>*Method</i>	*Automated		*Automated		*Manual		*Automated	
High Temp Alarm Setpoint = -20.0C	Alarms at _____C Temp flashes? Y N		Alarms at _____C Temp flashes? Y N		Alarms at _____C Temp flashes? Y N		Alarms at _____C Temp flashes? Y N	
Tech ID								
Date								
Reviewed by								
Date								

ANNUAL PM:

Probes (if necessary) Replace Battery	
Probe Bottles Dirty?	Y N
Probe Bottles Cleaned?	Y N
Tech ID	
Date	
Reviewed by	
Date	

AS NEEDED PM:

Task	Tech ID	Date	Tech ID	Date
Check & refill/replace probe bottle solution				
Clean Freezer Exterior				
Clean Freezer Interior				
Check the seal on Door Gasket				
Clean the Door Gaskets				
Calibrate the Chart Recorder				
Reviewed by:				