

Sensoscientific quick tips. 6.18.24

- Log in to website. There should be a quicklink on the lab shared drive to do this.

GO TO: <https://cloud.sensoscientific.com/Account/Login.aspx>

USERNAME: your e-mail address

PASSWORD: GLHS1234

1. Once you are logged in click on CONFIGURE at the top of the page
 2. Click on USER in the orange box to the left of the screen
 3. Click on CHANGE PASSWORD in the gray box
 - a. Type in the old password GLHS1234
 - b. Type in your new password
 - c. Confirm your new password
 - d. Click on CHANGE PASSWORD
- Not everyone has been set up but if you feel you do need access, please let me know and I can add you to the system.
 - Not all lab temperature monitoring is hooked up to this. OP-LAB and MOB are the only off site that have temperature nodes currently. MEDC and CPH do still require manual temperatures to be taken which means thermometers need to be calibrated annually by the lab.
 - The lab is also helping to watch and monitor off site non-lab specimen fridges in Urology and Maria Stein. We also have the Occ Health drug screen fridge.
 - Alarms are set to go off when the temperature has been out for over 60 minutes. A phone call is made to the lab by an automated system.
 - Lab staff should verify errors online to see temp chart for item alarming.
 - Temp may be out for 3 reasons. 1) Range error 2) Connectivity error (Wi-Fi), 3) Battery error.
 - Techs should put reason and action information online to document what was done and if contents needed to be moved or if monitoring until further action is needed.
 - If offsite after hours, please comment and leave Tom or someone a message to call the office in the morning. Facilities may need to check on it.
 - Charts have been updated in a new format to record the status of fridge or recording of temperature. The only way to see trends is to go online and review the chart of the item in question.