

Procedure Change/Revision (IMM 36-A)

Procedure: Tecan Freedom EVO 100

Change/Revision:

Maintenance Revision: To be added to the *weekly maintenance* schedule.

1. Cleaning of the waster chute
 - a. Remove waste chute from disposal apparatus.
 - b. Soak in hot soapy water for a few hours or even overnight.
 - c. Use a toothbrush or Q-tip applicator stick to clean precipitated serum from the chute.
 - d. Dry thoroughly with a sports towel.

Date of Revision: 09/07/12

Director Signature/Date:

 9/4/2012

Technologist Signature/Date:

Refer to MTS.