




# ODH New BSL III Decontamination (Steriplex)

By Eric St Germain

# Steriplex

- ODH new disinfectant by sBioMed
- Steriplex is a 2 part system
- It must be activated before use and once activated is best if used within 60 days
- Inactivated- each part is stable until the expiration date on the container

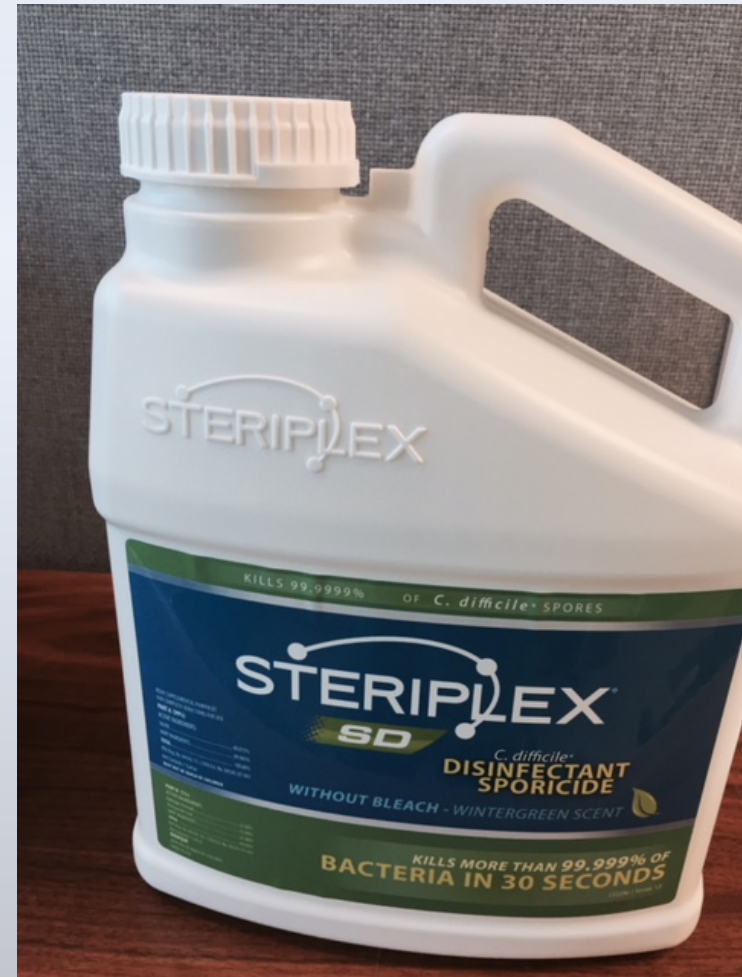

# Steriplex Activation

## Procedure

1. Remove the cap.
2. Push firmly on the number 1 location on the cap.
3. Place cap back on container and shake vigorously for 15 seconds.
4. Remove the cap.
5. Peel the film off lid. Steriplex is now ready to be used.
6. Write the date activated on the container and expiration date (60 days from activation)
7. Prepare 2 cans of wipes



# Wipe Preparation

## Procedure

Items needed: 2 new rolls of wipes, 2 labeled BioMed wipe containers, 1 L graduated cylinder and activated Steriplex.

1. Remove lid of the Wipe containers
2. Place new wipe roll into each container.

Follow wipe guidance diagram



# Wipe Preparation

## How to use Sanotex® Environmental Surface Wipes

A



Insert a dry roll of Sanotex® wipes into the canister.

B



Carefully pour chemical solution over center of wipe roll in canister.  
Allow five minutes before use to assure proper saturation of wipes.

C



From the center of the roll, pull a wipe up and carefully thread through opening in canister lid (approximately one inch). Replace lid, screwing on tightly.

D



To dispense, open the hinged lid and pull the wipe through the opening at an angle. The spring-loaded lid will re-close automatically.

# Wipe Preparation

## Procedure


3. Add 1 L of Activated Steriplex, pour into the center of the wipe roll.
4. From the center of the roll pull up on a wipe and place in the lid.
5. Screw on the lid and let it sit for 5 minutes before use.



# Steriplex Activation

## Procedure

6. On the back of the wipe containers write the date the Steriplex was activated. And the date of expiration (60 days).

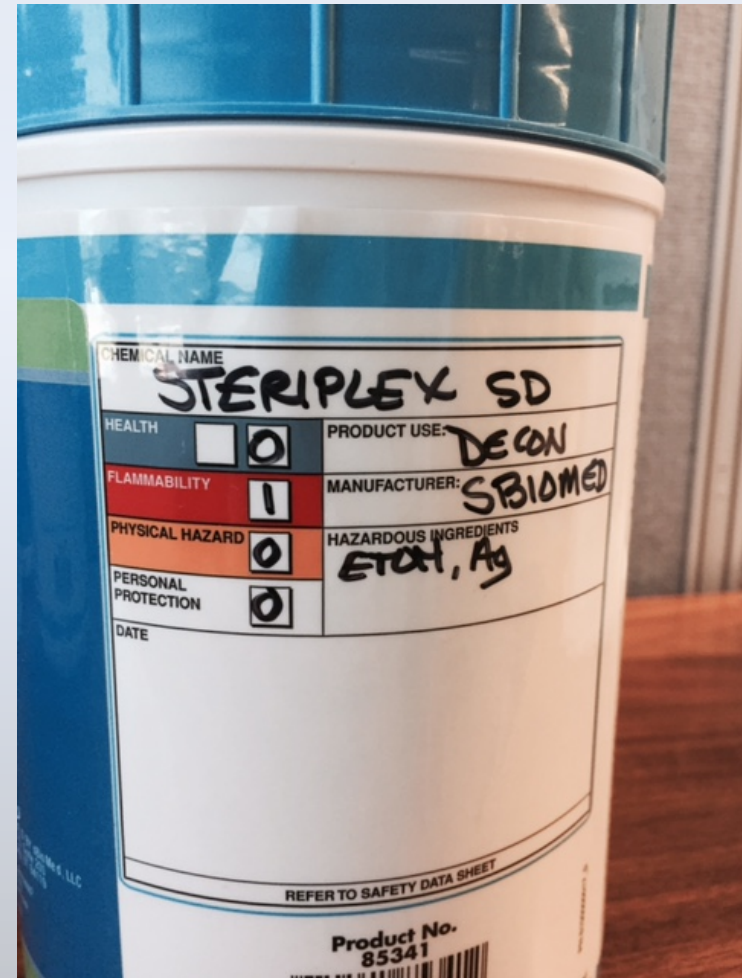

# Steriplex Activation

## Procedure

6. On the back of the wipe containers write the date the Steriplex was activated. And the date of expiration (60 days).

Please see the lead BT Trainer if you have any questions.

