


# Change Summary: Updated Shipper for BMS Cell Therapy Drug Products

BMS Confidential Information. © 2026 Bristol-Myers Squibb Company.  
03/2026 HE-US-2600245

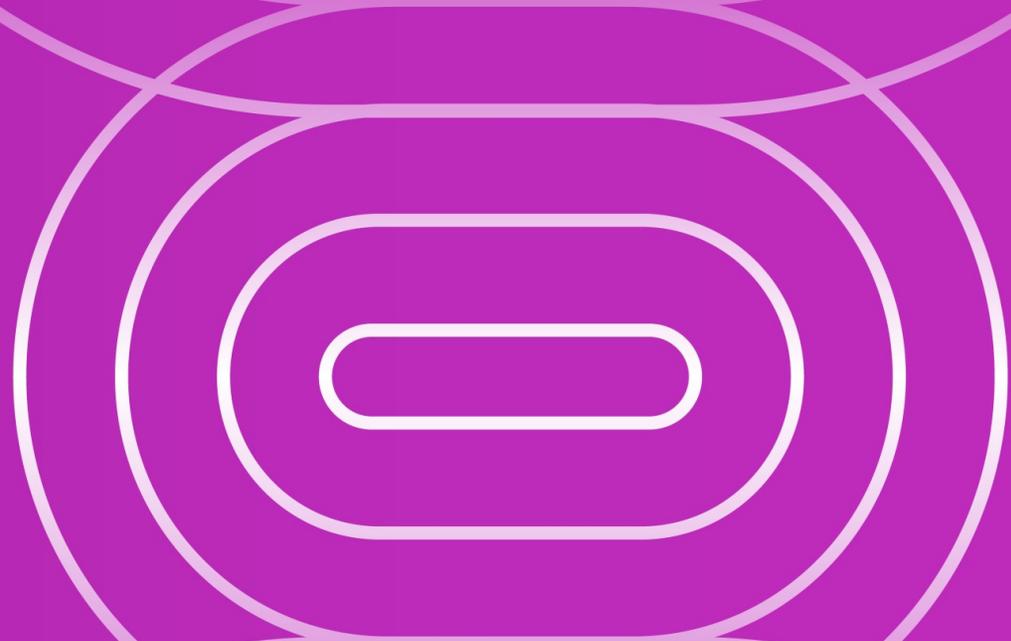

# Disclaimers

- This presentation is intended for training on operational matters for implementing CAR T Cell therapy at Sites of Care.
- Please note the intent is to discuss operational considerations and not to promote any products.
- The content of this presentation and related conversations are subject to a confidentiality agreement and should not be disseminated to anyone not covered by the confidentiality agreement.

# Change Summary

- BMS is implementing the Cryoport HV3 General Purpose Shipping System, featuring a new exterior shipping case for cryogenic shippers
- There is no change to:
  - The interior shipper chamber which holds drug product
  - Thermal performance or validation of the cryogenic shipper
  - Data logger or temperature monitoring process
  - Product handling procedures
- Change will be implemented starting 1 Apr 2026

## Current shipper:

- Cylindrical outer shell with rounded top
- Shipping container requires separate dolly or mechanism for transport
- Hanging document pouches inside
- Paper instructions included
- Cannot be stacked or palletized



## Updated Shipper (as of April 1, 2026)

- Cube-shaped outer shell with built-in dolly system
- Integrated document compartment (no hanging pouches)
- Designated exterior label panel
- Stackable design for transport
- Same interior chamber with 8-day validated hold time at cryogenic temperatures as current shipper

# Exterior Shipper Specifications

Product Specifications	Updated shipper as of April 1, 2026- HV3 Shipping System	Current shipper- HV2 Shipping System
Image		
Temperature Range	-150C or colder	-150C or colder
Dynamic Hold Time	Up to 8 days	Up to 8 days
Exterior Dimensions	22" L x 22" W x 28" H 558.8 mm x 558.8 mm x 711.2 mm	22" D x 28" H 558.8 mm x 711.2 mm
Interior Chamber Dimensions	8.5" (215.9 mm) diameter 12" (304.8 mm) height Max commodity height 11.5"	8.5" (215.9 mm) diameter 12" (304.8 mm) height Max commodity height 11.5"
Dry Weight	Up to 73 lb / 33 kg	Up to 57lb / 26 kg
Charged Weight	Up to 92 lb / 42 kg	Up to 77 lb / 35 kg

# Exterior Images



# Exterior Images

Waybills & shipping labels

Dolly handle



Shipper expiry label panel



Zip tie and latch closure

# Opening the shipper

1 Cut the zip-tie on each latch



2 Depress the button on the underside of the latch and lift to release




Push up button  
on the underside  
of the latch



# Document layout change - inside

## Current

### There Will Be Two Shipping Pouches Inside the Cryoshipper

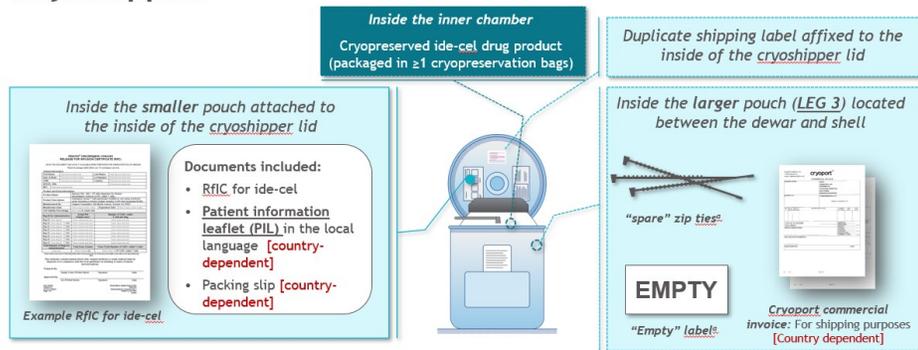

††treatment sites may decide whether to use the label, or zip ties. If the site does not need to use these items, the site may leave them untouched in the LEG 3 pouch.

PIL, patient information leaflet; RfIC, Release for Infusion Certificate.

References: 1. BMS internal files: 2022; 2. Abecma (idecabtagene vixtuocel) Release for Infusion Certificate. Varies from market to market.

Contact BMS at [PHONE NUMBER] with questions.

Bristol Myers Squibb | ide-cel Clinician Handling Guide

Subject to confidentiality agreement. Do not disseminate without prior authorization by BMS. | 28

## Future



2 plastic pouches included in the document compartment:

- Leg 2 will contain:
  - Release for Infusion Certificate
  - Patient information leaflet [country dependent]
  - Packing slip [country dependent]
  - **Secondary BMS shipper expiry label**
- Leg 3 will contain:
  - Empty label
  - Spare zip ties
  - Commercial invoice [country dependent]

# Accessing the Document Compartment




**WARNING:** The lid opens and rests at a 90-degree angle. Close it carefully to avoid slamming, as pushing it will cause it to shut abruptly.

Document compartment is clearly visible and accessible when the exterior lid is open

# Using the Built in Dolly to Move the Shipper

- 1 Slide lever to unlock handle and pull handle up with your other hand.

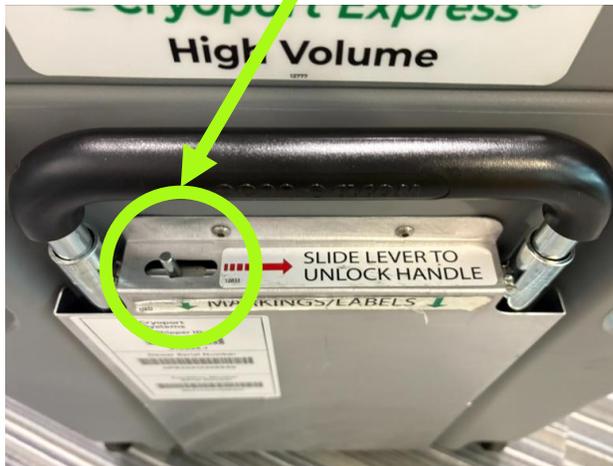

- 2 Lift handle until you hear an audible click. Handle can extend to two different heights. Tilt the shipper and use wheels to pull like a suitcase.



# Label Placement

Label requirements may vary by country





UN3245 sticker

Shipper expiry label form with fields for Shipper name, Full Delivery address for lab, and other details.

Shipper expiry label



Do Not X-Ray sticker [country dependent]

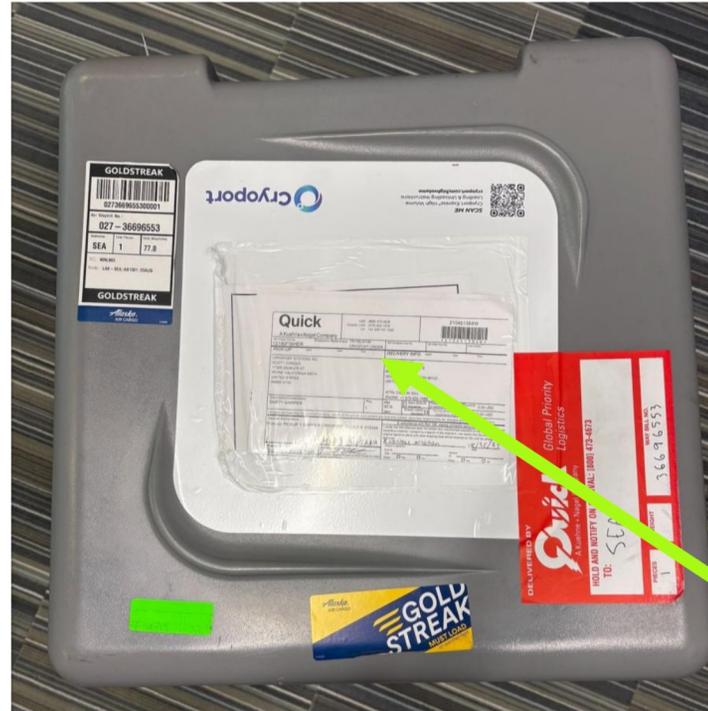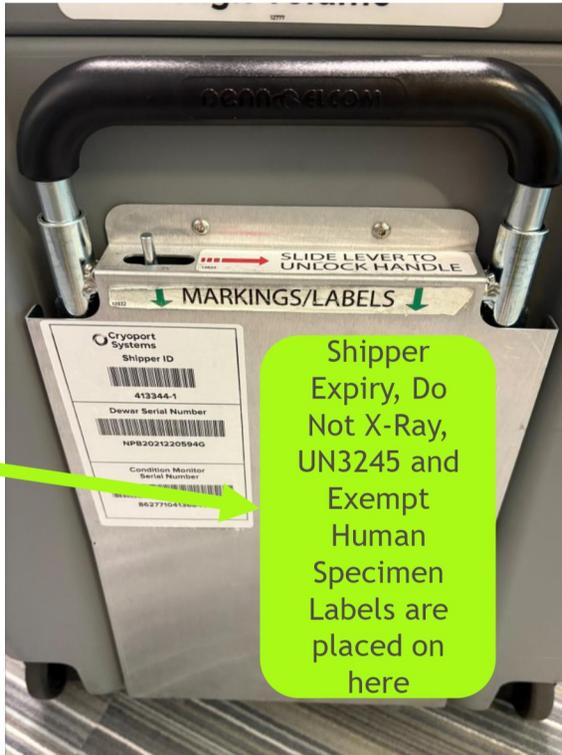

Waybill

Waybill and return label [country dependent] are placed here

# Return Shipper Instructions: Secure the latch with zipties



Thread the zip tie through the hole and close it around the latch




# Follow BMS Handling Instructions



ALWAYS FOLLOW  
BMS INSTRUCTIONS  
FOR CELL THERAPY  
DRUG PRODUCT AND  
SHIPPER HANDLING.

DISREGARD  
CRYOPORT.COM  
INSTRUCTIONS  
LINKED TO AT THIS  
QR CODE.

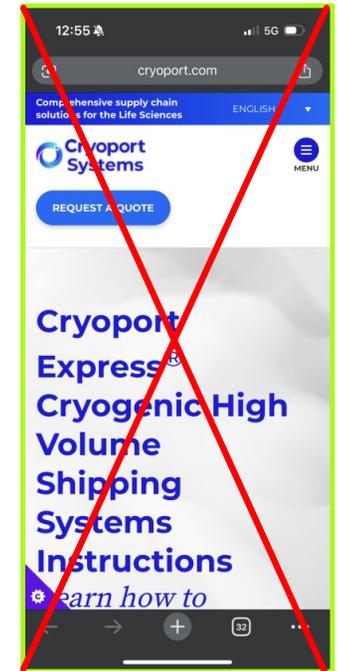

# Contact BMS CT360 Scheduling for any questions

[Scheduling@celltherapy360.com](mailto:Scheduling@celltherapy360.com)